


Computer Controlled Multiple Compressor Refrigeration Control, with SCADA







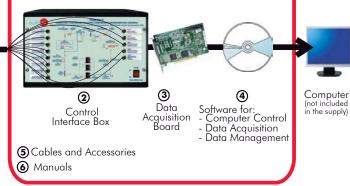

EDIBON SCADA System

1 Unit: THALAC/1. Multiple Compressor Refrigeration Control

*Minimum supply always includes: 1 + 2 + 3 + 4 + 5 + 6 (Computer not included in the supply)

Key features:

- Advanced Real-Time SCADA.
- Open Control + Multicontrol + Real-Time Control.
- Specialized EDIBON Control Software based on Labview.
- National Instruments Data Acquisition board (250 KS/s, kilo samples per second).
- Calibration exercises, which are included, teach the user how to calibrate a sensor and the importance of checking the accuracy of the sensors before taking measurements.
- ▶ Projector and/or electronic whiteboard compatibility allows the unit to be explained and demonstrated to an entire class at one time.
- Capable of doing applied research, real industrial simulation, training courses, etc.
- Remote operation and control by the user and remote control for EDIBON technical support, are always included.
- Totally safe, utilizing 4 safety systems (Mechanical, Electrical, Electronic & Software).
- Designed and manufactured under several quality standards.
- ▶ Optional CAL software helps the user perform calculations and comprehend the results.
- This unit has been designed for future expansion and integration. A common expansion is the EDIBON Scada-Net (ESN) System which enables multiple students to simultaneously operate many units in a network.

For more information about Key Features, click here:







Certificates ISO 14000 and ECO-Management and Audit Scheme (environmental management)

www.edibon.com

Products
Products range
Units

3-Thermodynamics
8 Thermotechnics

OPEN CONTROL

MULTICONTROL

REAL TIME CONTROL







INTRODUCTION -

In traditional refrigeration units the cooling efficiency is controlled by including more condensers to reduce the temperature, meaning a higher efficiency in the evaporator. The problem of this type of control is that the compressor is always working; therefore, the electricity consumption is the same although at that moment the needs do not require so much consumption.

Other way of controlling the cooling efficiency is including more compressors, so that when only one compressor is working, the electricity consumption decreases considerably.

The objective of the Multiple Compressor Refrigeration Control (THALAC/1) unit is to introduce the student into the complex world of heat pumps systems, as well as studying and determining the characteristic operation parameters of the unit in function of the environmental requirements (humidity, heat, temperature and cooling, etc.).

GENERAL DESCRIPTION

The THALAC/1 has been designed to allow the students to understand the behaviour and characteristics of a heat pump (with air evaporator and water condenser) working in the steam compression cycle and with three compressors connected in parallel.

The refrigeration control unit with several compressors is basically composed of a refrigeration circuit. There is a water condenser, an air evaporator and three compressors located along it.

The low-grade heat source will be the air in the atmosphere. The heat generated by this unit is in form of hot water.

The work introduced is in form of electric energy supplied to the hermetic compressors.

This Computer Controlled Unit is supplied with the EDIBON Computer Control System (SCADA), and includes: The unit itself + a Control Interface Box + a Data Acquisition Board + Computer Control, Data Acquisition and Data Management Software Packages, for controlling the process and all parameters involved in the process.

PROCESS DIAGRAM AND UNIT ELEMENTS ALLOCATION 4 actuators and 12 sensors **OPEN CONTROL** controlled from any computer, MULTICONTROL and working simultaneously REAL TIME CONTROL FLOW SENSOR TEMPERATURE SENSOR **TEMPERATURE SENSOR** MANOMETER MANOMETER Sensor de Caudal Sensor de temperatura Manómetro Sensor de temperatura Manómetro FILTER **EXPANSION VALVE** LIQUID RECEIVER Válvula de Expansión Acumulador **Filtro** AEAI-1 AEA-1 AIR EVAPORATOR WATER CONDENSER Condensador de Agua ST-6, **Evaporador de Aire** ST-8 PRESSURE SENSOR Sensor de Presión ACO-3 COMPRESSOR Compresor SP-1 ST-4 TEMPERATURE SENSOR **TEMPERATURE SENSOR** PRESSURE Sensor de temperatura Sensor de temperatura SENSOR ACO-2 Sensor de Presión COMPRESSOR M-4 MANOMETER Compresor MANOMETER Manómetro Manómetro LIQUID SEPARATOR Separador de ACO-1 AP-1 líquido **SW-1 HIGH PRESSURE** COMPRESSOR WATTMETER Compresor Sensor de Potencia

Page 2 www.edibon.com

COMPLETE TECHNICAL SPECIFICATIONS (for main items)

With this unit there are several options and possibilities:

- Main items: 1, 2, 3, 4, 5 and 6.
- Optional items: 7, 8, 9, 10, 11 and 12.

Let us describe first the main items (1 to 6):

① THALAC/1. Unit:

Bench-top unit.

Anodized aluminum structure and panels in painted steel.

Main metallic elements in stainless steel.

Diagram in the front panel with similar distribution to the elements in the real unit.

3 Cooling compressors, computer controlled:

Technical data of each compressor:

Power: 1/5CV, 1.9A.

Fridge efficiency: 460W to 0°C. Cylinder volume: 7.55 cm³.

This compound system is controlled so that individual compressor can be

switched depending on the performance.

Water condenser:

Concentric pipes type. Capacity: 1580W. Consume DT=8K:170 I./h.

Coolant accumulation tank, capacity: 1.50 dm³.

Cooling filter.

Expansion valve, capillar type.

Air evaporator, computer controlled. Air flow: 980 m³/h.

Tank of division of the cooling liquid.

2 Low pressure manometers.

2 High pressure manometers.

High pressure control:

Pressure switch: the pressure switch disconnect the electric power to the compressor when the pressure in the circuit is higher than the fixed pressure.

Type: Spring, with automatic restart.

Switch off pressure: 14 Bar.

Restart pressure: 9.5 Bar.

8 Temperature sensors "J" type:

Refrigerant temperature in the compressor outlet.

Refrigerant temperature in the condenser outlet.

Refrigerant temperature in the evaporator inlet.

Refrigerant temperature in the compressor inlet.

Water temperature in the condenser inlet.

Water temperature in the condenser outlet.

Environmental temperature.

Air temperature in the evaporator outlet.

2 Flow sensors:

Cooling flow sensor, magnetic type, range: 5 to 60 l./h.

Water flow sensor (condenser), turbine type, range: 0.2 to 6.5 l./min.

2 Pressure sensors:

High pressure sensor, range: 0 to 25 bar.

Low pressure sensor, range: 0 to 10 bar.

Power measurement form the computer (PC).

Enthalpy diagram of the refrigerant R134a.

The complete unit includes as well:

Advanced Real-Time SCADA.

Open Control + Multicontrol + Real-Time Control.

Specialized EDIBON Control Software based on Labview.

National Instruments Data Acquisition board (250 KS/s, kilo samples per second).

Calibration exercises, which are included, teach the user how to calibrate a sensor and the importance of checking the accuracy of the sensors before taking measurements.

Projector and/or electronic whiteboard compatibility allows the unit to be explained and demonstrated to an entire class at one time.

Capable of doing applied research, real industrial simulation, training courses, etc.

Remote operation and control by the user and remote control for EDIBON technical support, are always included.

Totally safe, utilizing 4 safety systems (Mechanical, Electrical, Electronic & Software).

Designed and manufactured under several quality standards.

Optional CAL software helps the user perform calculations and comprehend the results.

This unit has been designed for future expansion and integration. A common expansion is the EDIBON Scada-Net (ESN) System which enables multiple students to simultaneously operate many units in a network.



THALAC/1. Unit

Page 3

② THALAC/1/CIB. Control Interface Box:

The Control Interface Box is part of the SCADA system.

<u>Control interface box with process diagram in the front panel</u> and with the same distribution that the different elements located in the unit, for an easy understanding by the student.

All sensors, with their respective signals, are properly manipulated from -10V. to +10V. computer output. Sensors connectors in the interface have different pines numbers (from 2 to 16), to avoid connection errors.

Single cable between the control interface box and computer.

The unit control elements are permanently computer controlled, without necessity of changes or connections during the whole process test procedure.

Simultaneous visualization in the computer of all parameters involved in the process.

Calibration of all sensors involved in the process.

Real time curves representation about system responses.

Storage of all the process data and results in a file.

Graphic representation, in real time, of all the process/system responses.

All the actuators' values can be changed at any time from the keyboard allowing the analysis about curves and responses of the whole process.

All the actuators and sensors values and their responses are displayed on only one screen in the computer.

Shield and filtered signals to avoid external interferences.

Real time computer control with flexibility of modifications from the computer keyboard of the parameters, at any moment during the process.

Real time computer control for pumps, compressors, resistances, control valves, etc.

Real time computer control for parameters involved in the process simultaneously.

Open control allowing modifications, at any moment and in real time, of parameters involved in the process simultaneously.

Three safety levels, one mechanical in the unit, another electronic in the control interface and the third one in the control software.

3 DAB. Data Acquisition Board:

The Data Acquisition board is part of the SCADA system.

PCI Express Data acquisition board (National Instruments) to be placed in a computer slot.

Bus PCI Express.

Analog input:

Number of **channels = 16** single-ended or 8 differential. **Resolution = 16 bits**, 1 in 65536.

Sampling rate up to: 250 KS/s (kilo samples per second).

Input range (V) = ± 10 V. Data transfers = DMA, interrupts, programmed I/0. DMA channels = 6.

Analog output:

Number of channels=2. Resolution=16 bits, 1 in 65536. Maximum output rate up to: 900 KS/s.

Output range(V) = ± 10 V. Data transfers = DMA, interrupts, programmed I/0.

Digital Input/Output:

Number of channels=24 inputs/outputs. D0 or DI Sample Clock frequency: 0 to 100 MHz.

Timing: Number of Counter/timers=4. Resolution: Counter/timers: 32 bits.

@ THALAC/1/CCSOF. Computer Control + Data Acquisition + Data Management Software:

The three softwares are part of the SCADA system.

Compatible with actual Windows operating systems.

Graphic and intuitive simulation of the process in screen.

Compatible with the industry standards.

Registration and visualization of all process variables in an automatic and simultaneous way.

Flexible, open and multicontrol software, developed with actual windows graphic systems, acting simultaneously on all process parameters.

Management, processing, comparison and storage of data.

Sampling velocity up to 250 KS/s (kilo samples per second).

Calibration system for the sensors involved in the process.

It allows the registration of the alarms state and the graphic representation in real time.

Comparative analysis of the obtained data, after the process and modification of the conditions during the process.

Open software, allowing the teacher to modify texts, instructions. Teacher's and student's passwords to facilitate the teacher's control on the student, and allowing the access to different work levels.

This unit allows the 30 students of the classroom to visualize simultaneously all the results and the manipulation of the unit, during the process, by using a projector or an electronic whiteboard.

⑤ Cables and Accessories, for normal operation.

Manuals: This unit is supplied with 8 manuals: Required Services, Assembly and Installation, Interface and Control Software, Starting-up, Safety, Maintenance, Calibration & Practices Manuals.

* References 1 to 6 are the main items: THALAC/1 + THALAC/1/CIB + DAB + THALAC/1/CCSOF + Cables and Accessories + Manuals are included in the minimum supply for enabling normal and full operation.



THALAC/1/CIB



SUCIONAL SENSOR AND SE

THALAC/1/CCSOF

EXERCISES AND PRACTICAL POSSIBILITIES TO BE DONE WITH MAIN ITEMS

1.- Combined operation of compressors:

Power measurement.

Comparison of the energy for operating individual compressor and multiple compressors.

- 2.- Cyclic process on the p-h state diagram.
- 3.- Determination of the inlet power, heat produced and performance coefficient. Air as heat source.
- Preparation of performance curves of the unit with different inlet and outlet temperatures. Air as heat source.
- 5.- Lay out of the steam compression cycle in a diagram P-H and comparison with the ideal cycle. Air as heat source.
- 6.- Effect of refrigerant supercooling.
- 7.- Effect of the airflow rate on the condenser performance.
- 8.- Preparation of the performance curves of the unit based on the properties of the refrigerant and at different condensation and evaporation temperatures. Air as heat source.
- 9.- Energy balances.

Other possible practices:

- 10.- Temperature sensors calibration.
- 11.- Flow sensors calibration.
- 12.- Pressure sensors calibration.

Other possibilities to be done with this Unit:

13.-Many students view results simultaneously.

To view all results in real time in the classroom by means of a projector or an electronic whiteboard.

14.-Open Control, Multicontrol and Real Time Control.

This unit allows intrinsically and/or extrinsically to change the span, gains; proportional, integral, derivate parameters; etc, in real time.

- 15.-The Computer Control System with SCADA allows a real industrial simulation.
- 16.-This unit is totally safe as uses mechanical, electrical and electronic, and software safety devices.
- 17.-This unit can be used for doing applied research.
- 18.- This unit can be used for giving training courses to Industries even to other Technical Education Institutions.
- 19.-Control of the THALAC/1 unit process through the control interface box without the computer.
- Visualization of all the sensors values used in the THALAC/1 unit process.
- By using PLC-Pl additional 19 more exercises can be done.
- Several other exercises can be done and designed by the user.

REQUIRED SERVICES

- Electrical supply: 220V., 1-phase + neutral + ground, 50Hz ; or 110V., 1-phase + neutral + ground, 60Hz.; and 1CV max.
- Water supply and drainage.
- Computer (PC).

DIMENSIONS & WEIGHTS

THALAC/1:

Unit: -Dimensions: 1000 x 600 x 700 mm. approx.

(39.37 x 23.62 x 27.55 inches approx.

-Weight: 100 Kg. approx.

(220 pounds approx.)

Control-Interface Box: -Dimensions: 490 x 330 x 310 mm. approx.

(19.29 x 12.99 x 12.2 inches approx.)

-Weight: 10 Kg. approx.

(22 pounds approx.)


RECOMMENDED ACCESSORIES

For refilling R134a refrigerant and maintenance, we recommend:

- T/KIT1. Maintenance Kit, containing: vacuum pump, hoses and manometers.
- T/KIT2. Maintenance Kit, containing: leakage detector.
- R134a refrigerant (to be acquired by the customer locally).

Page 5 www.edibon.com


SCADA Main screen

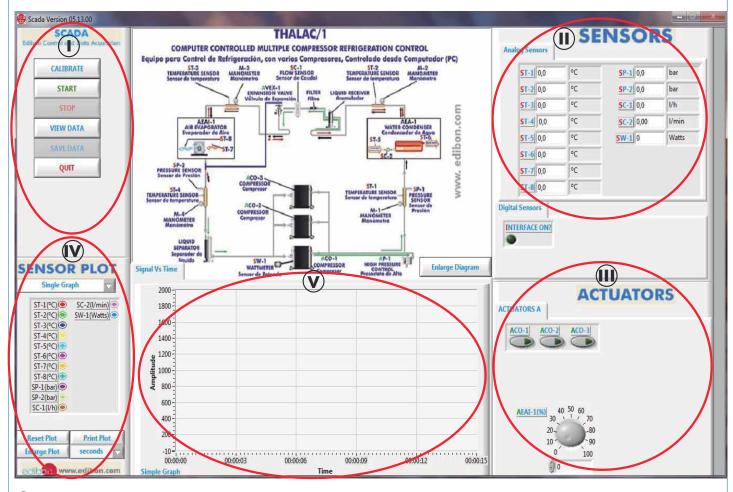

- Main software operation possibilities.
- (II) Sensors displays, real time values, and extra output parameters. Sensors: ST=Temperature sensor. SP=Pressure sensor. SC=Flow sensor. SW= Wattmeter.
- (ii) Actuators controls. Actuators: ACO= Compressor. AEAI= Air evaporator.
- (V) Channel selection and other plot parameters.
- Real time graphics displays.

Software for Sensors Calibration


Example of screen

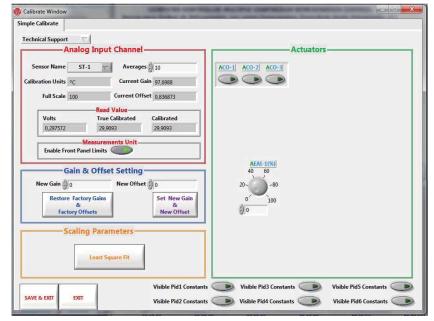

By using a free of charge code, the teacher and the students can calibrate the unit. The teacher can recover his/her own calibration by using the EDIBON code that we give free of charge.


Page 6 www.edibon.com

Using one compressor:

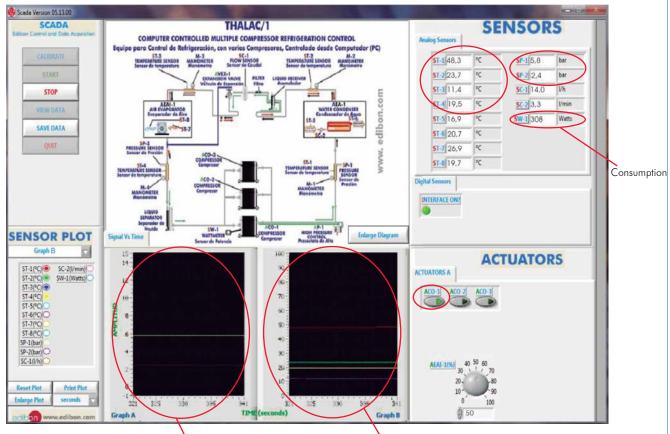


High and low pressures (SP-1, SP-2)

Refrigeration circuit temperatures (ST-1, ST-2, ST-3, ST-4)

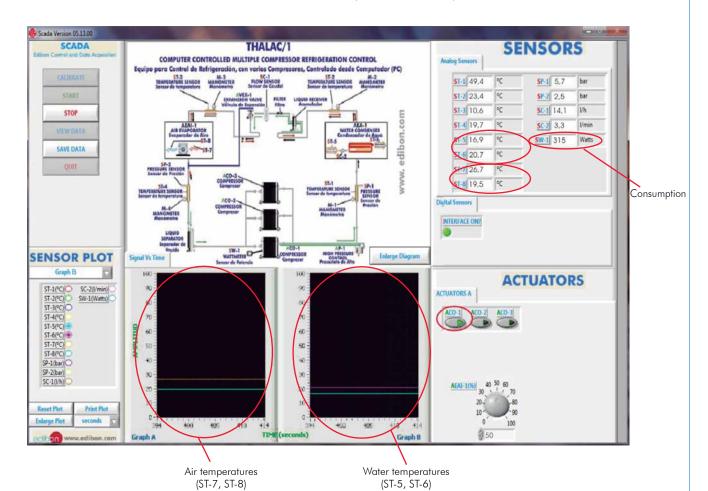


Continue...

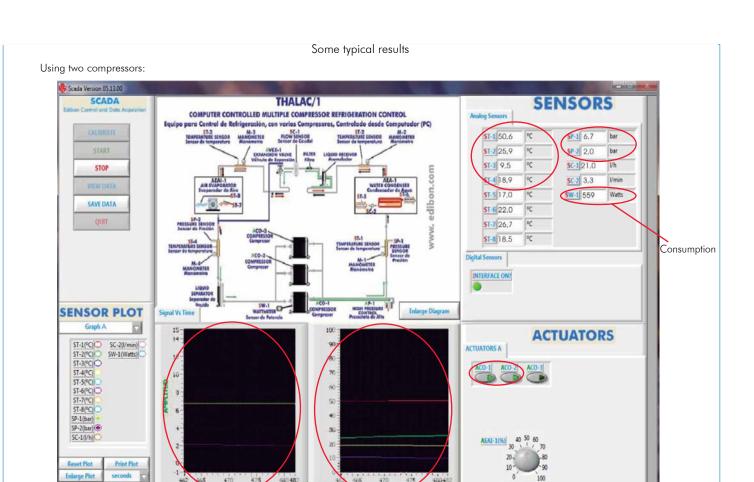

TIM

Pressures

(SP-1, SP-2)

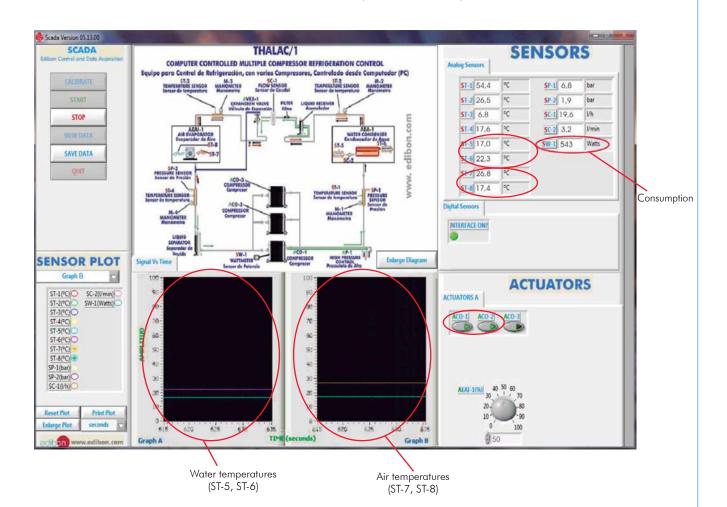

Page 8

Refrigeration circuit temperatures

(ST-1, ST-2, ST-3, ST-4)

Continue...

Some typical results

Using three compressors:


Reset Plot


Print Plot

Water temperatures (ST-5, ST-6)

The THALAC/1 unit allows the activation and operation of 3 compressors connected in parallel. It allows to study the system with one compressor and to compare its efficiency with that obtained when working with 3 compressors. The software allows to regulate (0-100%) the fan that works as evaporator.

When the three compressors are used a greater coolant flow rate, as well as a smaller temperature difference between ST-1 and ST-2 and a greater temperature difference between ST-3 and ST-4, can be observed.

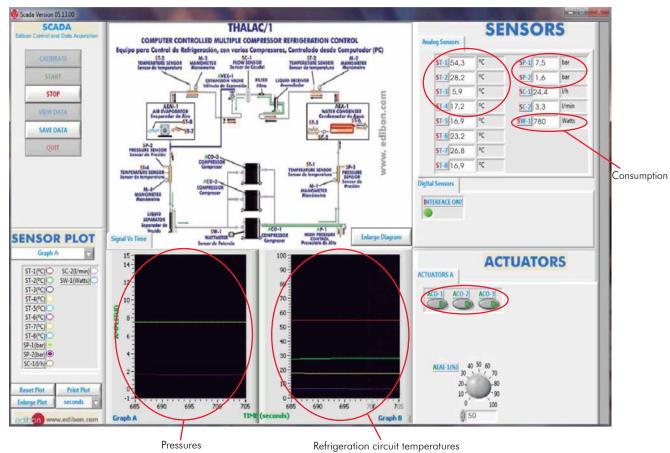

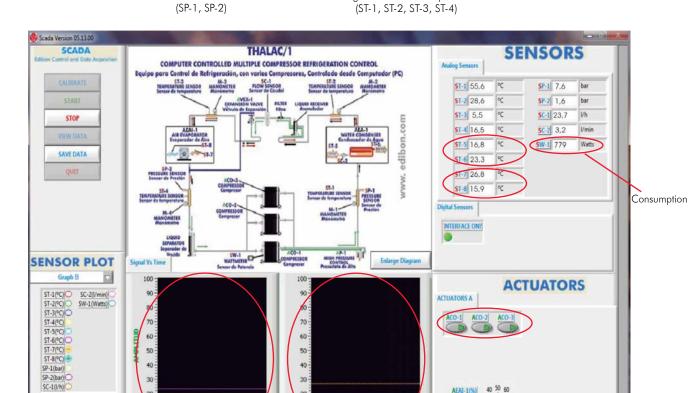

760

Air temperatures (ST-7, ST-8)

Page 9 www.edibon.com

COMPLETE TECHNICAL SPECIFICATIONS (for optional items)

Additionally to the main items (1 to 6) described, we can offer, as optional, other items from 7 to 12.

All these items try to give more possibilities for:

- a) Industrial configuration. (PLC)
- b) Technical and Vocational Education configuration. (CAI and FSS)
- c) Higher Education and/or Technical and Vocational Education configuration. (CAL)
- d) Multipost Expansions options. (Mini ESN and ESN)

a) Industrial configuration

PLC. Industrial Control using PLC (it includes PLC-PI Module plus PLC-SOF Control Software):

-PLC-PI. PLC Module:

Metallic box.

Circuit diagram in the module front panel.

Front panel:

Digital inputs(X) and Digital outputs (Y) block:

16 Digital inputs, activated by switches and 16 LEDs for confirmation (red).

14 Digital outputs (through SCSI connector) with 14 LEDs for message (green).

Analog inputs block:

16 Analog inputs (-10 V. to + 10 V.) (through SCSI connector).

Analog outputs block:

4 Analog outputs (-10 V. to + 10 V.) (through SCSI connector).

Touch screen:

High visibility and multiple functions. Display of a highly visible status. Recipe function. Bar graph function. Flow display function. Alarm list. Multi language function. True type fonts.

Back panel:

Power supply connector. Fuse 2A. RS-232 connector to PC. USB 2.0 connector to PC.

Inside:

Power supply outputs: 24 Vdc, 12 Vdc, -12 Vdc, 12 Vdc variable.

Panasonic PLC:

High-speed scan of 0.32 μ sec. for a basic instruction.

Program capacity of 32 Ksteps, with a sufficient comment area.

Power supply input (100 to 240 V AC).

DC input: 16 (24 V DC).

Relay output: 14.

High-speed counter.

Multi-point PID control.

Digital inputs/outputs and analog inputs/outputs Panasonic modules.


Communication RS232 wire to computer (PC).

Dimensions: 490 x 330 x 310 mm. approx. (19.29 x 12.99 x 12.20 inches approx.). Weight: 30 Kg. approx. (66 pounds approx.).

-THALAC/1/PLC-SOF. PLC Control Software:

For this particular unit, always included with PLC supply.

The software has been designed using Labview and it follows the unit operation procedure and linked with the Control Interface Box used in the Computer Controlled Multiple Compressor Refrigeration Control (THALAC/1).

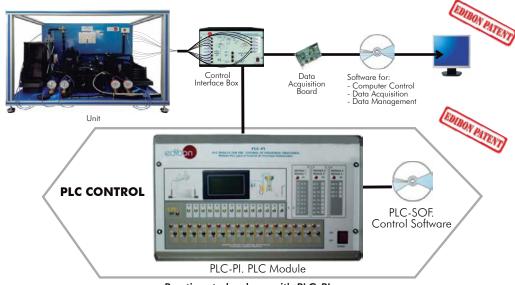

Practices to be done with PLC-PI:

- Control of the THALAC/1 unit process through the control interface box without the computer.
- 2.- Visualization of all the sensors values used in the THALAC/1 unit process.
- 3.- Calibration of all sensors included in the THALAC/1 unit process.
- 4.- Hand on of all the actuators involved in the THALAC/1 unit process.
- 5.- Realization of different experiments, in automatic way, without having in front the unit. (This experiment can be decided previously).
- 6.- Simulation of outside actions, in the cases hardware elements do not exist. (Example: test of complementary tanks, complementary industrial environment to the process to be studied, etc).
- 7.- PLC hardware general use and manipulation.
- 8.- PLC process application for THALAC/1 unit.
- 9.- PLC structure.

- 10.- PLC inputs and outputs configuration.
- 11.- PLC configuration possibilities.
- 12.- PLC programming languages.
- 13.- PLC different programming standard languages.
- 14.- New configuration and development of new process.
- 15.- Hand on an established process.
- 16.- To visualize and see the results and to make comparisons with the $\mathsf{THALAC}/1$ unit process.
- 17.- Possibility of creating new process in relation with the THALAC/1 unit.
- 18.- PLC Programming exercises.
- 19.- Own PLC applications in accordance with teacher and student requirements.

b) Technical and Vocational Education configuration

® THALAC/1/CAI. Computer Aided Instruction Software System.

This complete software package includes two Softwares: the INS/SOF. Classroom Management Software (Instructor Software) and the THALAC/1/SOF. Computer Aided Instruction Software (Student Software).

This software is optional and can be used additionally to items (1 to 6).

This complete software package consists on an Instructor Software (INS/SOF) totally integrated with the Student Software (THALAC/1/SOF). Both are interconnected so that the teacher knows at any moment what is the theoretical and practical knowledge of the students.

- INS/SOF. Classroom Management Software (Instructor Software): The Instructor can:

Organize Students by Classes and Groups.

Create easily new entries or delete them.

Create data bases with student information.

Analyze results and make statistical comparisons.

Generate and print reports.

Detect student's progress and difficulties.

...and many other facilities.

-THALAC/1/SOF. Computer Aided Instruction Software (Student Software): It explains how to use the unit, run the experiments and what to do at any moment.

This Software contains:

Theory.

Exercises

Guided Practices.

Exams.

For more information see CAI catalogue. Click on the following link:

www.edibon.com/products/catalogues/en/CAI.pdf

CAL CAL INSTRUCTOR SOFTWARE



Instructor Software

Student Software



Example of some screens

THALAC/1/FSS. Faults Simulation System.

Faults Simulation System (FSS) is a Software package that simulates several faults in any EDIBON Computer Controlled Unit. It is useful for Technical and Vocational level.

The "FAULTS" mode consists on causing several faults in the unit normal operation. The student must find them and solve them.

There are several kinds of faults that can be grouped in the following sections:

Faults affecting the sensors measurement:

- An incorrect calibration is applied to them.
- Non-linearity.

Faults affecting the actuators:

- Actuators channels interchange at any time during the program
- Response reduction of an actuator.

Faults in the controls execution:

- Inversion of the performance in ON/OFF controls.
- Reduction or increase of the calculated total response.
- The action of some controls is annulled.

On/off faults:

- Several on/off faults can be included.

For more information see FSS catalogue. Click on the following link: www.edibon.com/products/catalogues/en/FSS.pdf

Page 11 www.edibon.com

c) Higher Education and/or Technical and Vocational Education configuration

① THALAC/1/CAL. Computer Aided Learning Software (Results Calculation and Analysis).

This Computer Aided Learning Software (CAL) is a Windows based software, simple and very easy to use, specifically developed by EDIBON. It is very useful for Higher Education level.

CAL is a class assistant that helps in doing the necessary calculations to extract the right conclusions from data obtained during the experimental

CAL computes the value of all the variables involved and performs the calculations.

It allows to plot and print the results. Within the plotting options, any variable can be represented against any other.

Different plotting displays.


It has a wide range of information, such as constant values, unit conversion factors and integral and derivative tables.

For more information see CAL catalogue. Click on the following link:

www.edibon.com/products/catalogues/en/CAL.pdf



Information of constant values, unit conversion factors and integral and derivative tables

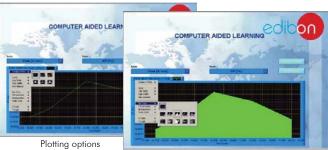

d) Multipost Expansions options

① Mini ESN. EDIBON Mini Scada-Net System.

Mini ESN. EDIBON Mini Scada-Net System allows up to 30 students to work with a Teaching Unit in any laboratory, simultaneously. It is useful for both, Higher Education and/or Technical and Vocational Education.

The Mini ESN system consists on the adaptation of any EDIBON computer controlled unit with SCADA integrated in a local network.

This system allows to view/control the unit remotely, from any computer integrated in the local net (in the classroom), through the main computer connected to the unit. Then, the number of possible users who can work with the same unit is higher than in an usual way of working (usually only one).

Main characteristics:

- It allows up to 30 students to work simultaneously with the EDIBON Computer Controlled Unit with SCADA, connected in a local net.
- Open Control + Multicontrol + Real Time Control + Multi Student Post.
- Instructor controls and explains to all students at the same time.
- Any user/student can work doing "real time" control/multicontrol and visualisation.
- Instructor can see in the computer what any user/student is doing in the unit.
- Continuous communication between the instructor and all the users/ students connected.

Main advantages:

- It allows an easier and quicker understanding.
- This system allows you can save time and cost.
- Future expansions with more EDIBON Units.

For more information see Mini ESN catalogue. Click on the following link: www.edibon.com/products/catalogues/en/Mini-ESN.pdf

Mini ESN. **EDIBON Mini Scada-Net System** 1 UNIT = up to 30 STUDENTS can work simultaneously Multiple Compressor Refrigeration Control (THALAC/1) Computer Control Software: Computer Control+Data Control Interface Acquisition+Data Management nstructor Central Mini Scada-Net Software 30 Student Posts LOCAL NET **OPEN CONTROL** Note: The Mini ESN MULTICONTROL system can be used with any EDIBON REAL TIME CONTROL computer controlled MULTI STUDENT POST

@ESN. EDIBON Scada-Net System.

This unit can be integrated, in the future, into a Complete Laboratory with many Units and many Students.

For more information see **ESN** catalogue. Click on the following link:

www.edibon.com/products/catalogues/en/units/thermodynamicsthermotechnics/esn-thermodynamics/ESN-THERMODYNAMICS.pdf

Page 12 www.edibon.com

ORDER INFORMATION

Main items (always included in the supply)

Minimum supply always includes:

- ① Unit: THALAC/1. Multiple Compressor Refrigeration Control.
- ② THALAC/1/CIB. Control Interface Box.
- ③ DAB. Data Acquisition Board.
- 4 THALAC/1/CCSOF. Computer Control + Data Acquisition + Data Management Software.
- ⑤ Cables and Accessories, for normal operation.
- Manuals.
- * <u>IMPORTANT:</u> Under <u>THALAC/1</u> we always supply all the elements for immediate running as 1, 2, 3, 4, 5 and 6.

Optional items (supplied under specific order)

a) Industrial configuration

- PLC. Industrial Control using PLC (it includes PLC-PI Module plus PLC-SOF Control Software):
 - PCL-PI. PLC Module.
 - THALAC/1/PLC-SOF. PLC Control Software.

b) Technical and Vocational configuration

- ❸ THALAC/1/CAI. Computer Aided Instruction Software System.
- c) <u>Higher Education and/or Technical and Vocational Education configuration</u>
- THALAC/1/CAL. Computer Aided Learning Software (Results Calculation and Analysis).

d) Multipost Expansions options

- 10 Mini ESN. EDIBON Mini Scada-Net System.
- 19 ESN. EDIBON Scada-Net System.

Page 13 www.edibon.com

TENDER SPECIFICATIONS (for main items)

①THALAC/1. Unit:

Bench-top unit.

Anodized aluminum structure and panels in painted steel.

Main metallic elements in stainless steel

Diagram in the front panel with similar distribution to the elements in the real unit.

3 Cooling compressors, computer controlled:

Technical data of each compressor:

Power: 1/5CV, 1.9A.

Fridge efficiency: 460W to 0°C.

Cylinder volume: 7.55 cm³

This compound system is controlled so that individual compressor can be switched depending on the performance.

Water condenser:

Concentric pipes type. Capacity: 1580W. Consume DT=8K:170 I./h.

Coolant accumulation tank, capacity: 1.50 dm³.

Cooling filter.

Expansion valve, capillar type.

Air evaporator, computer controlled. Air flow: 980 m³/h.

Tank of division of the cooling liquid.

2 Low pressure manometers.

2 High pressure manometers

High pressure control:

Pressure switch: the pressure switch disconnect the electric power to the compressor when the pressure in the circuit is higher than the fixed pressure .

Type: Spring, with automatic restart.

Switch off pressure: 14 Bar

Restart pressure: 9.5 Bar.

8 Temperature sensors "J" type:

Refrigerant temperature in the compressor outlet.

Refrigerant temperature in the condenser outlet.

Refrigerant temperature in the evaporator inlet.

Refrigerant temperature in the compressor inlet.

Water temperature in the condenser inlet.

Water temperature in the condenser outlet.

Environmental temperature.

Air temperature in the evaporator outlet.

2 Flow sensors:

Cooling flow sensor, magnetic type, range: 5 to 60 l./h.

Water flow sensor (condenser), turbine type, range: 0.2 to 6.5 l./min.

2 Pressure sensors:

High pressure sensor, range: 0 to 25 bar.

Low pressure sensor, range: 0 to 10 bar.

Power measurement form the computer (PC).

Enthalpy diagram of the refrigerant R134a.

The complete unit includes as well:

Advanced Real-Time SCADA.

Open Control + Multicontrol + Real-Time Control.

Specialized EDIBON Control Software based on Labview

National Instruments Data Acquisition board (250 KS/s, kilo samples per second).

Calibration exercises, which are included, teach the user how to calibrate a sensor and the importance of checking the accuracy of the sensors before taking measurements.

Projector and/or electronic whiteboard compatibility allows the unit to be explained and demonstrated to an entire class at one time.

Capable of doing applied research, real industrial simulation, training courses, etc.

Remote operation and control by the user and remote control for EDIBON technical support, are always included

Totally safe, utilizing 4 safety systems (Mechanical, Electrical, Electronic & Software)

Designed and manufactured under several quality standards. Optional CAL software helps the user perform calculations and comprehend the results.

This unit has been designed for future expansion and integration. A common expansion is the EDIBON Scada-Net (ESN) System which enables multiple students to simultaneously operate many units in a network.

②THALAC/1/CIB. Control Interface Box:

The Control Interface Box is part of the SCADA system. Control interface box with process diagram in the front panel.

The unit control elements are permanently computer controlled.

Simultaneous visualization in the computer of all parameters involved in the process.

Calibration of all sensors involved in the process.

Real time curves representation about system responses.

All the actuators' values can be changed at any time from the keyboard allowing the analysis about curves and responses of the whole process.

Shield and filtered signals to avoid external interferences.

Real time computer control with flexibility of modifications from the computer keyboard of the parameters, at any moment during the process.

Real time computer control for parameters involved in the process simultaneously.

Open control allowing modifications, at any moment and in real time, of paraméters involved in the process simultaneously.

Three safety levels, one mechanical in the unit, another electronic in the control interface and the third one in the control software.

③ DAB. Data Acquisition Board:

The Data Acquisition board is part of the SCADA system.

PCI Express Data acquisition board (National Instruments) to be placed in a computer slot.

Analog input: Channels = 16 single-ended or 8 differential. Resolution = 16 bits, 1 in 65536. Sampling rate up to: 250 KS/s (kilo samples per second).

Analog output: Channels=2. Resolution=16 bits, 1 in 65536.

Digital Input/Output: Channels=24 inputs/outputs THALAC/1/CCSOF. Computer Control + Data Acquisition + Data Management Software:

The three softwares are part of the SCADA system.

Compatible with the industry standards.

Flexible, open and multicontrol software, developed with actual windows graphic systems, acting simultaneously on all process parameters.

Management, processing, comparison and storage of data. Sampling velocity up to 250 KS/s (kilo samples per second).

Calibration system for the sensors involved in the process.

It allows the registration of the alarms state and the graphic representation in real time.

Open software, allowing the teacher to modify texts, instructions. Teacher's and student's passwords to facilitate the teacher's control on the student, and allowing the access to different work levels.

This unit allows the 30 students of the classroom to visualize simultaneously all the results and the manipulation of the unit, during the process, by using a projector or an electronic whiteboard.

(5) Cables and Accessories, for normal operation.

6 Manuals: This unit is supplied with 8 manuals: Required Services, Assembly and Installation, Interface and Control Software, Starting-up, Safety, Maintenance, Calibration & Practices Manuals.

> Page 14 www.edibon.com

Exercises and Practical Possibilities to be done with Main Items

1.- Combined operation of compressors:

Power measurement.

Comparison of the energy for operating individual compressor and multiple compressors.

- 2.- Cyclic process on the p-h state diagram.
- 3.- Determination of the inlet power, heat produced and performance coefficient. Air as heat source.
- 4.- Preparation of performance curves of the unit with different inlet and outlet temperatures. Air as heat source.
- 5.- Lay out of the steam compression cycle in a diagram P-H and comparison with the ideal cycle. Air as heat source.
- 6.- Effect of refrigerant supercooling.
- 7.- Effect of the airflow rate on the condenser performance.
- 8.- Preparation of the performance curves of the unit based on the properties of the refrigerant and at different condensation and evaporation temperatures. Air as heat source.
- 9.- Energy balances.

Other possible practices:

- 10.- Temperature sensors calibration.
- 11.- Flow sensors calibration.
- 12.- Pressure sensors calibration.

Other possibilities to be done with this Unit:

13.- Many students view results simultaneously.

To view all results in real time in the classroom by means of a projector or an electronic whiteboard.

14.- Open Control, Multicontrol and Real Time Control.

This unit allows intrinsically and/or extrinsically to change the span, gains; proportional, integral, derivate parameters; etc, in real time.

- 15.- The Computer Control System with SCADA allows a real industrial simulation.
- 16.- This unit is totally safe as uses mechanical, electrical and electronic, and software safety devices.
- 17.- This unit can be used for doing applied research.
- 18.- This unit can be used for giving training courses to Industries even to other Technical Education Institutions.
- 19.- Control of the THALAC/1 unit process through the control interface box without the computer.
- 20.- Visualization of all the sensors values used in the THALAC/1 unit process.
- By using PLC-PI additional 19 more exercises can be done.
- Several other exercises can be done and designed by the user.

TENDER SPECIFICATIONS (for optional items)

a) Industrial configuration

PLC. Industrial Control using PLC (it includes PLC-PI Module plus PLC-SOF Control Software):

-PLC-PI. PLC Module:

Metallic box.

Circuit diagram in the module front panel.

Digital inputs(X) and Digital outputs (Y) block: 16 Digital inputs. 14 Digital outputs.

Analog inputs block: 16 Analog inputs.

Analog outputs block: 4 Analog outputs.

Touch screen.

Panasonic PLC:

 $High-speed\ scan\ of\ 0.32\ \mu sec.\ Program\ capacity\ of\ 32\ Ksteps.\ High-speed\ counter.\ Multi-point\ PID\ control.$

Digital inputs/outputs and analog inputs/outputs Panasonic modules.

-THALAC/1/PLC-SOF. PLC Control Software:

For this particular unit, always included with PLC supply.

Practices to be done with PLC-PI:

- 1.- Control of the THALAC/1 unit process through the control interface box without the computer.
- 2.- Visualization of all the sensors values used in the THALAC/1 unit process.
- 3.- Calibration of all sensors included in the THALAC/1 unit process.
- 4.- Hand on of all the actuators involved in the THALAC/1 unit process.
- 5. Realization of different experiments, in automatic way, without having in front the unit. (This experiment can be decided previously).
- 6.- Simulation of outside actions, in the cases hardware elements do not exist. (Example: test of complementary tanks, complementary industrial environment to the process to be studied, etc).
- 7.- PLC hardware general use and manipulation.
- 8.- PLC process application for THALAC/1 unit.
- 9.- PLC structure.
- 10.- PLC inputs and outputs configuration.
- 11.- PLC configuration possibilities.
- 12.- PLC programming languages.
- 13.- PLC different programming standard languages.
- 14.- New configuration and development of new process.
- 15.- Hand on an established process.
- 16.- To visualize and see the results and to make comparisons with the THALAC/1 unit process.
- 17.- Possibility of creating new process in relation with the THALAC/1 unit.
- 18.- PLC Programming exercises.
- $19.- Own PLC \ applications \ in \ accordance \ with \ teacher \ and \ student \ requirements.$

b) Technical and Vocational Education configuration

® THALAC/1/CAI. Computer Aided Instruction Software System.

This complete software package consists on an Instructor Software (INS/SOF) totally integrated with the Student Software (THALAC/1/SOF).

-INS/SOF. Classroom Management Software (Instructor Software):

The Instructor can:

Organize Students by Classes and Groups.

Create easily new entries or delete them.

Create data bases with student information.

Analyze results and make statistical comparisons.

Generate and print reports.

Detect student's progress and difficulties.

-THALAC/1/SOF. Computer Aided Instruction Software (Student Software):

It explains how to use the unit, run the experiments and what to do at any moment.

This Software contains:

Theory.

Exercises.

Guided Practices.

Exams.

THALAC/1/FSS. Faults Simulation System.

Faults Simulation System (FSS) is a Software package that simulates several faults in any EDIBON Computer Controlled Unit.

The "FAULTS" mode consists on causing several faults in the unit normal operation. The student must find them and solve them.

There are several kinds of faults that can be grouped in the following sections:

Faults affecting the sensors measurement:

- An incorrect calibration is applied to them.
- Non-linearity.

Faults affecting the actuators:

- Actuators channels interchange at any time during the program execution.
- Response reduction of an actuator.

Faults in the controls execution:

- Inversion of the performance in $\ensuremath{\mathsf{ON}}\xspace/\ensuremath{\mathsf{OFF}}\xspace$ controls.
- Reduction or increase of the calculated total response.
- The action of some controls is annulled.

On/off faults:

- Several on/off faults can be included.

Page 16 www.edibon.com

c) Higher Education and/or Technical and Vocational Education configuration

This Computer Aided Learning Software (CAL) is a Windows based software, simple and very easy to use.

CAL is a class assistant that helps in doing the necessary calculations to extract the right conclusions from data obtained during the experimental practices.

CAL computes the value of all the variables involved and performs the calculations.

It allows to plot and print the results. Within the plotting options, any variable can be represented against any other.

Different plotting displays.

It has a wide range of information, such as constant values, unit conversion factors and integral and derivative tables.

d) Multipost Expansions options

11) Mini ESN. EDIBON Mini Scada-Net System.

EDIBON Mini Scada-Net System allows up to 30 students to work with a Teaching Unit in any laboratory, simultaneously.

The Mini ESN system consists on the adaptation of any EDIBON computer controlled unit with SCADA integrated in a local network.

This system allows to view/control the unit remotely, from any computer integrated in the local net (in the classroom), through the main computer connected to the unit

Main characteristics:

- -It allows up to 30 students to work simultaneously with the EDIBON Computer Controlled Unit with SCADA, connected in a local net.
- -Open Control + Multicontrol + Real Time Control + Multi Student Post.
- -Instructor controls and explains to all students at the same time.
- -Any user/student can work doing "real time" control/multicontrol and visualisation.
- -Instructor can see in the computer what any user/student is doing in the unit.
- -Continuous communication between the instructor and all the users/students connected.

Main advantages:

- -It allows an easier a quicker understanding.
- -This system allows you can save time and cost.
- -Future expansions with more EDIBON Units.

The system basically will consist of:

This system is used with a Computer Controlled Unit.

- -Instructor's computer.
- -Students' computers.
- -Local Network.
- -Unit-Control Interface adaptation.
- -Unit Software adaptation.
- -Webcam.
- -Mini ESN Software to control the whole system.
- -Cables and accessories required for a normal operation.

*Specifications subject to change without previous notice, due to the convenience of improvements of the product.



C/Del Agua, 14. Polígono Industrial San José de Valderas. 28918 LEGANÉS. (Madrid). SPAIN.

Phone: 34-91-6199363 FAX: 34-91-6198647

E-mail: edibon@edibon.com WEB site: www.edibon.com

Issue: ED01/14 Date: April/2014

n	_	п	n	_	C	_	Ν	т	۸.	ΤI	١.	
ĸ	г	М	ĸ	г	٠.	г	$I \setminus$	ш	н	ш	·V	Е