



Computer Controlled Reversible Heat Pump + Air Conditioning + Refrigeration with 2 Condensers and 2 Evaporators (Water/Air), with SCADA

THIBAR22C

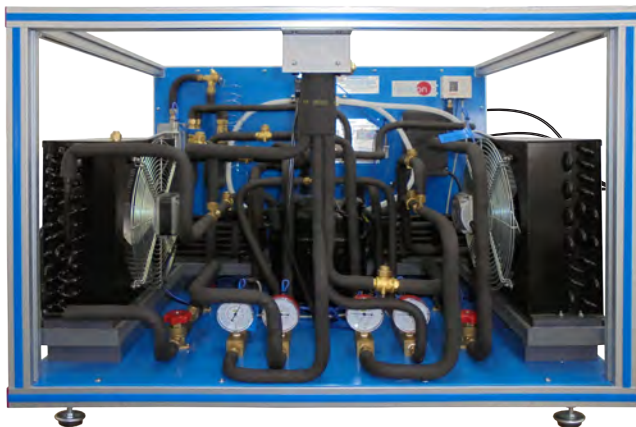

① Unit: THIBAR22C. Computer Controlled Reversible Heat Pump + Air Conditioning + Refrigeration with 2 Condensers and 2 Evaporators (Water/Air)

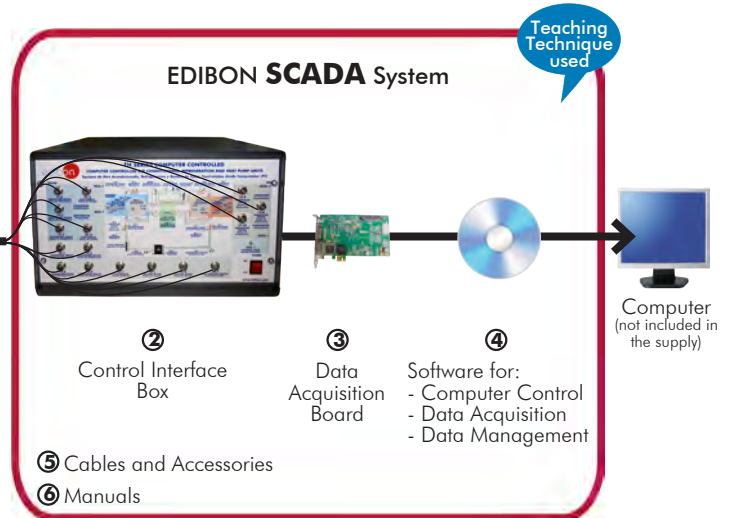

* Minimum supply always includes: 1 + 2 + 3 + 4 + 5 + 6
(Computer not included in the supply)

Key features:

- **Advanced Real-Time SCADA.**
- **Open Control + Multicontrol + Real-Time Control.**
- **Specialized EDIBON Control Software based on LabVIEW.**
- **National Instruments Data Acquisition board (250 KS/s, kilo samples per second).**
- **Calibration exercises, which are included, teach the user how to calibrate a sensor and the importance of checking the accuracy of the sensors before taking measurements.**
- **Projector and/or electronic whiteboard compatibility allows the unit to be explained and demonstrated to an entire class at one time.**
- **Capable of doing applied research, real industrial simulation, training courses, etc.**
- **Remote operation and control by the user and remote control for EDIBON technical support, are always included.**
- **Totally safe, utilizing 4 safety systems (Mechanical, Electrical, Electronic & Software).**
- **Designed and manufactured under several quality standards.**
- **Optional ICAI software to create, edit and carry out practical exercises, tests, exams, calculations, etc. Apart from monitoring user's knowledge and progress reached.**
- **This unit has been designed for future expansion and integration. A common expansion is the EDIBON Scada-Net (ESN) System which enables multiple students to simultaneously operate many units in a network.**

**OPEN CONTROL
+
MULTICONTROL
+
REAL TIME CONTROL**



www.edibon.com

PRODUCTS
↳ 9 - THERMODYNAMICS
& THERMOTECHNICS

For more information about Key Features, click here



ISO 9001: Quality Management (for Design, Manufacturing, Commercialization and After-sales service)



European Union Certificate (total safety)



Certificates ISO 14001 and ECO-Management and Audit Scheme (environmental management)



"Worlddidac Quality Charter" and Platinum Member of Worlddidac

INTRODUCTION

A heat pump is a device that provides thermal energy from a cold focus to hotter other one. For this, it is necessary a contribution of work, so in agreement with the 2^o law of the thermodynamics, the heat is transferred in a spontaneous way from a hot focus to cold other one, until their temperatures are equal.

The Computer Controlled Rversible Heat Pump + Air Conditioning + Refrigeration with 2 Condensers and 2 Evaporators (Water/Air), "THIBAR22C", is an unit designed to study the thermodynamic processes that happen in a circuit that it can work as heat pump, air conditioning and refrigeration and for which a coolant circulates.

GENERAL DESCRIPTION

The Computer Controlled Rversible Heat Pump + Air Conditioning + Refrigeration with 2 Condensers and 2 Evaporators (Water/Air), "THIBAR22C", has as aim introduce the student to the study of the heat pumps, air conditioning and refrigeration, as well as the analysis and determination of the operation typical parameters of the unit depending on the two types of fluids used in the processes of evaporation and condensation (air and water).

This unit can have different applications, depending on the type of cold focus or hot focus used in the processes of evaporation and condensation. This unit consists of the following stages:

Compression:

This stage begins when the coolant enters to the compressor. This coolant is compressed, increasing its pressure and temperature. To measure these variables the unit includes a pressure sensor, a manometer and a temperature sensor.



THIBAR22C detail

Condensation:

The coolant has two possibilities: divert the coolant through the air condenser, or through the water condenser. The coolant transfers its heat to the water (or to the air) that flows for the condenser. At the end of this stage, the pressure and the temperature of the coolant are measured by means of a manometer and a temperature sensor.

Expansion:

The coolant circulates through an accumulator and a filter, to retain particles of condensate, and a flow sensor. Next it circulates through the valve of expansion, which causes a fall of pressure and temperature of the coolant. At the end of this stage, the pressure and the temperature of the coolant are measured by means of a manometer and temperature sensor.

Evaporation:

The coolant has two possibilities: divert the coolant through the air evaporator, or through the water evaporator. The coolant absorbs the heat of the water (or the air) that flows for the evaporator. At the end of this stage, the pressure and the temperature of the coolant are measured by means of a pressure sensor, a manometer and a temperature sensor. Finally, the coolant circulates through a liquid separator to retain liquid particles before going on to the compressor.

The condensers and evaporators have different sensors to the measure of the most important parameters (temperatures and flows). In addition, the unit includes a high pressure control to avoid an excess of pressure in the unit and a four-way valve to change the direction of the coolant and to achieve the reversibility of the cycle.


The four-way valve (or cycle inversion valve) allows to obtain different combinations of Heat Pump, Air Conditioning and Refrigeration in a only unit. So that using the cycle inversion valve, we have nine units in one, six of them as Heat Pump + Air Conditioning + Refrigeration and three of them Heat Pump + Air Conditioning.

This Computer Controlled Unit is supplied with the EDIBON Computer Control System (SCADA), and includes: The unit itself + a Control Interface Box + a Data Acquisition Board + Computer Control, Data Acquisition and Data Management Software Packages, for controlling the process and all parameters involved in the process.

PROCESS DIAGRAM AND UNIT ELEMENTS ALLOCATION

4 actuators and 18 sensors controlled from any computer, and working simultaneously

OPEN CONTROL
+
MULTICONTROL
+
REAL TIME CONTROL

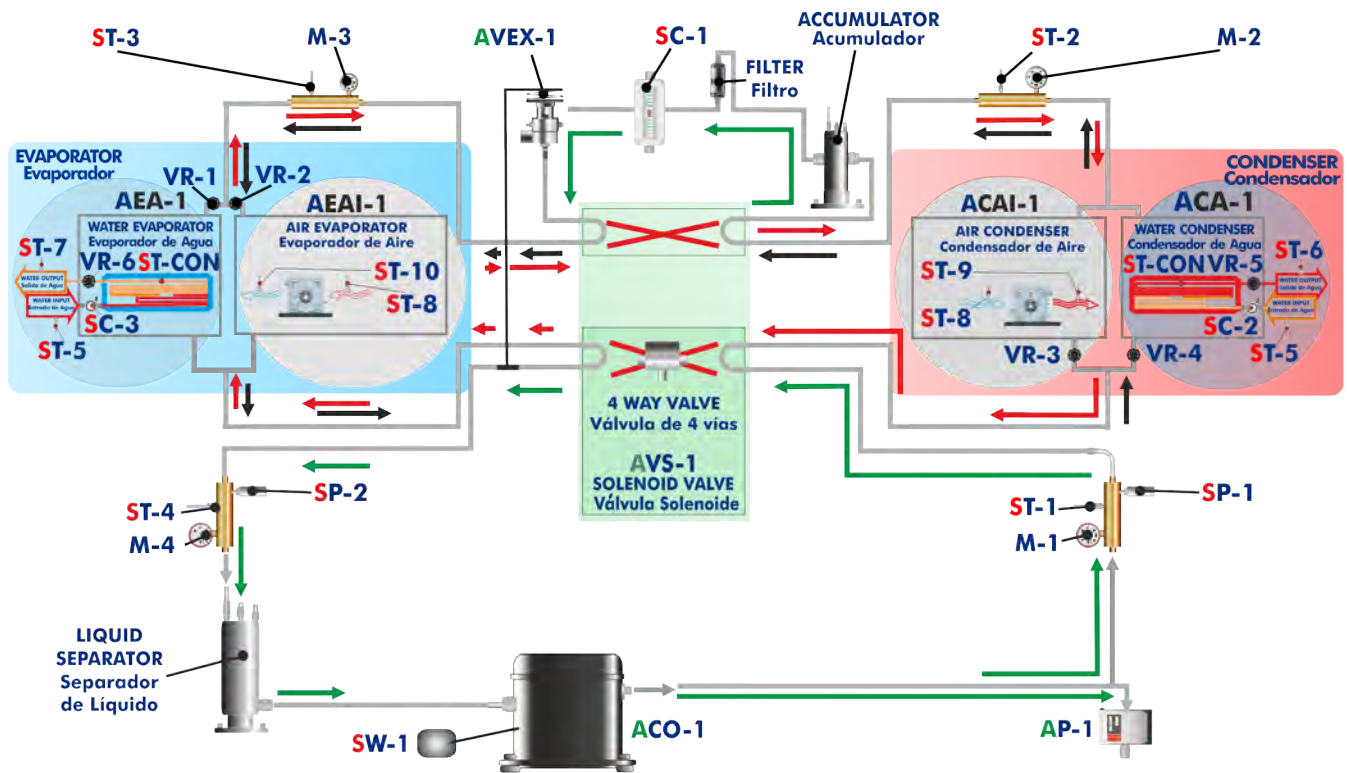

ST- TEMPERATURE SENSOR Sensor de Temperatura	SC- FLOW SENSOR Sensor de Caudal	M- MANOMETER Manómetro	SW- WATTMETER Sensor de Potencia	AP- HIGH PRESSURE CONTROL Presostato de Alta	VR- REGULATION VALVE Válvula de Regulación
ACO- COMPRESSOR Compresor	AVEX- EXPANSION VALVE Válvula de Expansión	ST-CON CONTROL TEMPERATURE SENSOR Sensor de Control de Temperatura	SP- PRESSURE SENSOR Sensor de Presión		

COMPLETE TECHNICAL SPECIFICATIONS (for main items)

With this unit there are several options and possibilities:

- Main items: 1, 2, 3, 4, 5 and 6.
- Optional items: 7, 8, 9, 10 and 11.

Let us describe first the main items (1 to 6):

① THIBAR22C. Unit:

Bench-top unit.

Anodized aluminum frame and panels made of painted steel.

Main metallic elements made of stainless steel.

Diagram in the front panel with distribution of the elements similar to the real one.

Hermetic compressor, power: 3/8 HP, computer controlled.

Water condenser, capacity: 1580 W.

Air condenser, computer controlled.

High pressure control.

Storage vessel to prevent a wrong operation of the compressor, capacity: 0.5 l.

Filter for the working fluid.

Expansion valve.

Water evaporator, capacity: 1580 W.

Air evaporator, computer controlled.

Liquid separator to prevent liquid from entering the compressor, capacity: 0.73 l.

Four manometers, situated at important points of the unit.

Ten temperature sensors (four sensors measure the cooling temperature, three sensors measure the water temperature, three sensors measure the air temperature):

- “J” type temperature sensor (compressor outlet).
- “J” type temperature sensor (condenser outlet/evaporator inlet).
- “J” type temperature sensor (evaporator inlet/condenser outlet).
- “J” type temperature sensor (compressor inlet).
- “J” type temperature sensor (water inlet).
- “J” type temperature sensor (condenser water outlet).
- “J” type temperature sensor (evaporator water outlet).
- “J” type temperature sensor (air inlet).
- “J” type temperature sensor (condenser air outlet).
- “J” type temperature sensor (evaporator air outlet).

Three flow sensors:

Cooling flow sensor, range: 5 – 60 l/h.

Water flow sensor (water condenser), range: 0.25 – 6.5 l/min.

Water flow sensor (water evaporator), range: 0.25 – 6.5 l/min.

Two pressure sensors:

Cooling pressure sensor (compressor outlet), range: 0 – 25 bar.

Cooling pressure sensor (compressor inlet), range: 0 – 10 bar.

Compressor wattmeter.

Four-way valve (solenoid valve) for reversibility the cycle.

Four valves to direct the coolant through the condensers and evaporators.

Enthalpy diagram of the R134a coolant.

The complete unit includes as well:

Advanced Real-Time SCADA.

Open Control + Multicontrol + Real-Time Control.

Specialized EDIBON Control Software based on LabVIEW.

National Instruments Data Acquisition board (250 KS/s, kilo samples per second).

Calibration exercises, which are included, teach the user how to calibrate a sensor and the importance of checking the accuracy of the sensors before taking measurements.

Projector and/or electronic whiteboard compatibility allows the unit to be explained and demonstrated to an entire class at one time.

Capable of doing applied research, real industrial simulation, training courses, etc.

Remote operation and control by the user and remote control for EDIBON technical support, are always included.

Totally safe, utilizing 4 safety systems (Mechanical, Electrical, Electronic & Software).

Designed and manufactured under several quality standards.

Optional ICAI software to create, edit and carry out practical exercises, tests, exams, calculations, etc. Apart from monitoring user's knowledge and progress reached.

This unit has been designed for future expansion and integration. A common expansion is the EDIBON Scada-Net (ESN) System which enables multiple students to simultaneously operate many units in a network.



Unit: THIBAR22C

② THIBAR22C/CIB. Control Interface Box:

The Control Interface Box is part of the SCADA system.

Control interface box with process diagram in the front panel and with the same distribution that the different elements located in the unit, for an easy understanding by the student.

All sensors, with their respective signals, are properly manipulated from -10V. to +10V. computer output. Sensors connectors in the interface have different pines numbers (from 2 to 16), to avoid connection errors.

Single cable between the control interface box and computer.

The unit control elements are permanently computer controlled, without necessity of changes or connections during the whole process test procedure.

Simultaneous visualization in the computer of all parameters involved in the process.

Calibration of all sensors involved in the process.

Real time curves representation about system responses.

Storage of all the process data and results in a file.

Graphic representation, in real time, of all the process/system responses.

All the actuators' values can be changed at any time from the keyboard allowing the analysis about curves and responses of the whole process.

All the actuators and sensors values and their responses are displayed on only one screen in the computer.

Shield and filtered signals to avoid external interferences.

Real time computer control with flexibility of modifications from the computer keyboard of the parameters, at any moment during the process.

Real time computer control for pumps, compressors, heating elements, control valves, etc.

Real time computer control for parameters involved in the process simultaneously.

Open control allowing modifications, at any moment and in real time, of parameters involved in the process simultaneously.

Three safety levels, one mechanical in the unit, another electronic in the control interface and the third one in the control software.



THIBAR22C/CIB

③ DAB. Data Acquisition Board:

The Data Acquisition board is part of the SCADA system.

PCI Express Data acquisition board (National Instruments) to be placed in a computer slot. Bus PCI Express.

Analog input:

Number of channels= 16 single-ended or 8 differential. Resolution=16 bits, 1 in 65536.

Sampling rate up to: 250 KS/s (kilo samples per second).

Input range (V)= ±10 V. Data transfers=DMA, interrupts, programmed I/O. DMA channels=6.

Analog output:

Number of channels=2. Resolution=16 bits, 1 in 65536.

Maximum output rate up to: 900 KS/s.

Output range (V)= ±10 V. Data transfers=DMA, interrupts, programmed I/O.

Digital Input/Output:

Number of channels=24 inputs/outputs. D0 or DI Sample Clock frequency: 0 to 100 MHz.

Timing: Number of Counter/timers=4. Resolution: Counter/timers: 32 bits.

The Data Acquisition board model may change at any moment, providing the same or better features than those required for the unit.



DAB

④ THIBAR22C/CCSOF. Computer Control + Data Acquisition + Data Management Software:

The three softwares are part of the SCADA system.

Compatible with actual Windows operating systems. Graphic and intuitive simulation of the process in screen. **Compatible with the industry standards.**

Registration and visualization of all process variables in an automatic and simultaneous way.

Flexible, open and multicontrol software, developed with actual windows graphic systems, acting simultaneously on all process parameters.

Management, processing, comparison and storage of data.

Sampling velocity up to 250 KS/s (kilo samples per second).

Calibration system for the sensors involved in the process.

It allows the registration of the alarms state and the graphic representation in real time.

Comparative analysis of the obtained data, after the process and modification of the conditions during the process.

Open software, allowing the teacher to modify texts, instructions. Teacher's and student's passwords to facilitate the teacher's control on the student, and allowing the access to different work levels.

This unit allows the 30 students of the classroom to visualize simultaneously all the results and the manipulation of the unit, during the process, by using a projector or an electronic whiteboard.



THIBAR22C/CCSOF

⑤ Cables and Accessories, for normal operation.

⑥ Manuals:

This unit is **supplied with 8 manuals**: Required Services, Assembly and Installation, Interface and Control Software, Starting-up, Safety, Maintenance, Calibration & Practices Manuals.

*References 1 to 6 are the main items: THIBAR22C + THIBAR22C/CIB + DAB + THIBAR22C/CCSOF + Cables and Accessories + Manuals are included in the minimum supply for enabling normal and full operation.

EXERCISES AND PRACTICAL POSSIBILITIES TO BE DONE WITH THE MAIN ITEMS

- | | |
|---|--|
| <ol style="list-style-type: none"> 1.- Determination of COP (coefficient of performance) of a heat pump. Water as heat source. (Water-water heat pump). 2.- Determination of COP (coefficient of performance) of a heat pump. Air as heat source. (Water-air heat pump). 3.- Determination of COP (coefficient of performance) of a heat pump. Air as heat source. (Air-air heat pump). 4.- Determination of COP (coefficient of performance) of a heat pump. Water as heat source. (Air-water heat pump). 5.- Representation of performance curves of the heat pump with different inlet and outlet temperatures. Water as heat source. (Water-water heat pump). 6.- Representation of performance curves of the heat pump with different inlet and outlet temperatures. Air as heat source. (Water-air heat pump). 7.- Representation of performance curves of the heat pump with different inlet and outlet temperatures. Water as heat source. (Air-water heat pump). 8.- Representation of performance curves of the heat pump with different inlet and outlet temperatures. Air as heat source. (Air-air heat pump). 9.- Lay out of the refrigeration compression cycle in a diagram P-H and comparison with the ideal cycle. Water as heat source. (Water-water heat pump). 10.- Lay out of the refrigeration compression cycle in a diagram P-H and comparison with the ideal cycle. Air as heat source. (Water-air heat pump). 11.- Lay out of the refrigeration compression cycle in a diagram P-H and comparison with the ideal cycle. Water as heat source. (Air-water heat pump). 12.- Lay out of the refrigeration compression cycle in a diagram P-H and comparison with the ideal cycle. Air as heat source. (Air-air heat pump). 13.- Representation of the performance curves of the heat pump based on the properties of the refrigerant and at different condensation and evaporation temperatures. Water as heat source. (Water-water heat pump). 14.- Representation of the performance curves of the heat pump based on the properties of the refrigerant and at different condensation and evaporation temperatures. Air as heat source. (Water-air heat pump). 15.- Representation of the performance curves of the heat pump based on the properties of the refrigerant and at different condensation and evaporation temperatures. Water as heat source. (Air-water heat pump). | <ol style="list-style-type: none"> 16.- Representation of the performance curves of the heat pump based on the properties of the refrigerant and at different condensation and evaporation temperatures. Air as heat source. (Air-air heat pump). 17.- Practices with cycle inversion. <p>Additional practical possibilities:</p> <ol style="list-style-type: none"> 18.- Sensors calibration. <p>Other possibilities to be done with this Unit:</p> <ol style="list-style-type: none"> 19.- Many students view results simultaneously.
To view all results in real time in the classroom by means of a projector or an electronic whiteboard. 20.- Open Control, Multicontrol and Real Time Control.
This unit allows intrinsically and/or extrinsically to change the span, gains; proportional, integral, derivative parameters; etc, in real time. 21.- The Computer Control System with SCADA allows a real industrial simulation. 22.- This unit is totally safe as uses mechanical, electrical/electronic, and software safety devices. 23.- This unit can be used for doing applied research. 24.- This unit can be used for giving training courses to Industries even to other Technical Education Institutions. 25.- Control of the THIBAR22C unit process through the control interface box without the computer. 26.- Visualization of all the sensors values used in the THIBAR22C unit process. <p>- By using PLC-PI additional 19 more exercises can be done.
- Several other exercises can be done and designed by the user.</p> |
|---|--|

REQUIRED SERVICES

- Electrical supply: single-phase 200 VAC – 240 VAC/50 Hz or 110 VAC – 127 VAC/60 Hz.
- Water supply and drain.
- Computer.

DIMENSIONS AND WEIGHTS

THIBAR22C:

Unit:

- Dimensions: 900 x 600 x 500 mm approx.
(35.43 x 23.62 x 19.68 inches approx.)
- Weight: 50 Kg approx.
(110 pounds approx.)

Control-Interface Box:

- Dimensions: 490 x 330 x 310 mm approx.
(19.29 x 12.99 x 12.20 inches approx.)
- Weight: 10 Kg approx.
(22 pounds approx.)

ADDITIONAL RECOMMENDED ELEMENTS (Not included)

For refilling R134a refrigerant and maintenance, we recommend:

- T/KIT1. Maintenance Kit, containing: vacuum pump, hoses and manometers.
- T/KIT2. Maintenance Kit, containing: leakage detector.
- R134a refrigerant (to be acquired by the customer locally).

SIMILAR UNITS AVAILABLE

Offered in this catalogue:

- THIBAR22C. Computer Controlled Rerversible Heat Pump + Air Conditioning + Refrigeration with 2 Condensers and 2 Evaporators (Water/Air).

Offered in other catalogue:

- THIBAR22B. Rerversible Heat Pump + Air Conditioning + Refrigeration with 2 Condensers and 2 Evaporators (Water/Air).
- THAR22C. Computer Controlled Heat Pump + Air Conditioning + Refrigeration with 2 Condensers and 2 Evaporators (Water/Air).
- THAR22B. Heat Pump + Air Conditioning + Refrigeration with 2 Condensers and 2 Evaporators (Water/Air).
- THIBAR44C. Computer Controlled Rerversible Heat Pump + Air Conditioning + Refrigeration with 4 Condensers and 4 Evaporators (Water/Air).
- THIBAR44B. Rerversible Heat Pump + Air Conditioning + Refrigeration with 4 Condensers and 4 Evaporators (Water/Air).

SOFTWARE MAIN SCREENS

SCADA Main screen

The SCADA main screen is divided into several functional areas:

- Top Left (I):** A control panel with buttons for CALIBRATE, START, STOP, VIEW DATA, SAVE DATA, and QUIT.
- Top Center:** A detailed schematic diagram of the THIBAR22C system, including components like CONDENSER, AIR CONDENSER, ACCUMULATOR, EVAPORATOR, and various valves and sensors.
- Top Right (II):** A 'SENSORS' display showing a grid of analog sensors (ST-1 to ST-10) and digital sensors (SP-1 to SP-2, SC-1 to SC-3, SW-1).
- Bottom Left (IV):** A 'SENSOR PLOT' window for 'SIGNAL VS TIME' with a list of sensors to plot and buttons for Reset Plot, Print Plot, and Enlarge Plot.
- Bottom Center (V):** A 'MOLLIER DIAGRAM' showing a graph of Amplitude vs. Time (seconds).
- Bottom Right (III):** An 'ACTUATORS' display showing controls for ACO-1 and AVS-1, with graphical representations of valves and compressors.

- (I) Main software operation possibilities.
- (II) Sensors displays, real time values, and extra output parameters. Sensors: ST= Temperature sensor. SP= Pressure sensor. SC= Flow sensor. SW= Wattmeter.
- (III) Actuators controls. Actuators: AVS= Solenoid valve (4-way valve). ACO= Compressor. AEAI= Air evaporator. ACAI= Air condenser.
- (IV) Channel selection and other plot parameters.
- (V) Real time graphics displays.

Software for Sensors Calibration

Example of screen

The 'Calibrate Window' interface is designed for sensor and actuator calibration. It features several key sections:

- Analog Input Channel:** Fields for Sensor Name (SC-2), Calibration Units (l/min), Full Scale (10), and Read Value (1.83321).
- Gain & Offset Setting:** Fields for New Gain (0) and New Offset (0), with buttons for 'Restore Factory Gains & Factory Offsets' and 'Set New Gain & New Offset'.
- Scaling Parameters:** A 'Least Square Fit' button.
- Actuators:** A 'Muticalibrate' section for ACO-1 and AVS-1, including graphical representations of the actuators.
- Sensors Table:** A detailed table for calibrating multiple sensors.


Sensors	Volts	Calibrated	Err(%)	Current Gain	Current Offset	Read	Set
SP-1	1,91531	9,75693	9,75693	5,06329	0,059157		
SP-2	1,9552	4,05233	4,05233	2,02154	0,09981		
SC-1	1,91282	19,8866	19,8866	10,4488	-0,1		
ST-1	1,94317	194,527	194,527	99,697	0,79894		
SW-1	1,91021	396,162	396,162	212,055	-8,9063		
ST-5	1,97892	194,648	194,648	97,3726	1,95597		
-	0	0	0	0	0		

Buttons at the bottom include 'Restore Factory Gains & Factory Offsets', 'Set New Gains & New Offsets', 'Take Measurement', 'Get New Gain & New Offset', and 'Taken Data'.


The teacher and the students can calibrate the unit with a password provided by EDIBON. The teacher can restore the factory calibration any time.

SOME REAL RESULTS OBTAINED FROM THIS UNIT

Real time representation of any measured variable. Evaporator and condenser water - water.

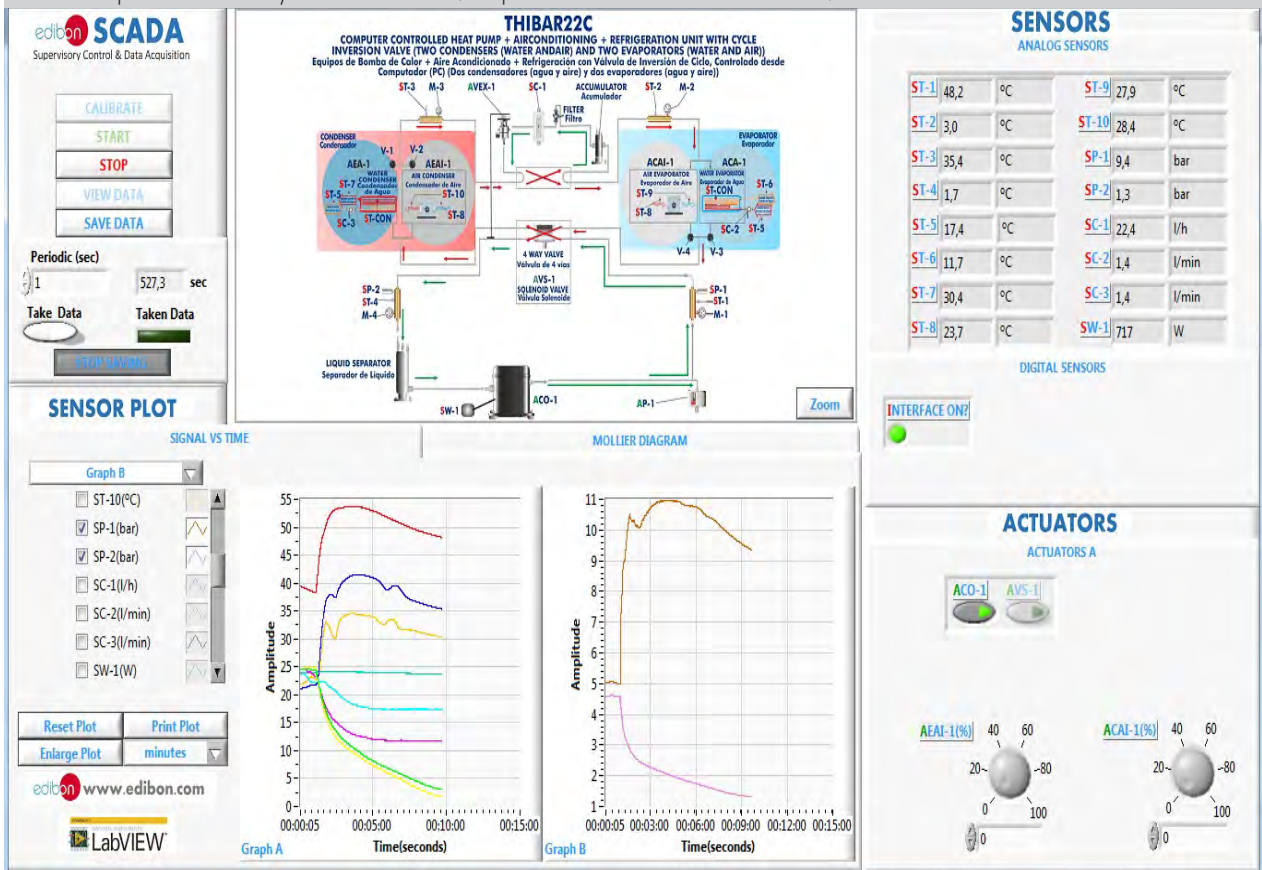

Real time representation of any measured variable. Evaporator and condenser air - air.

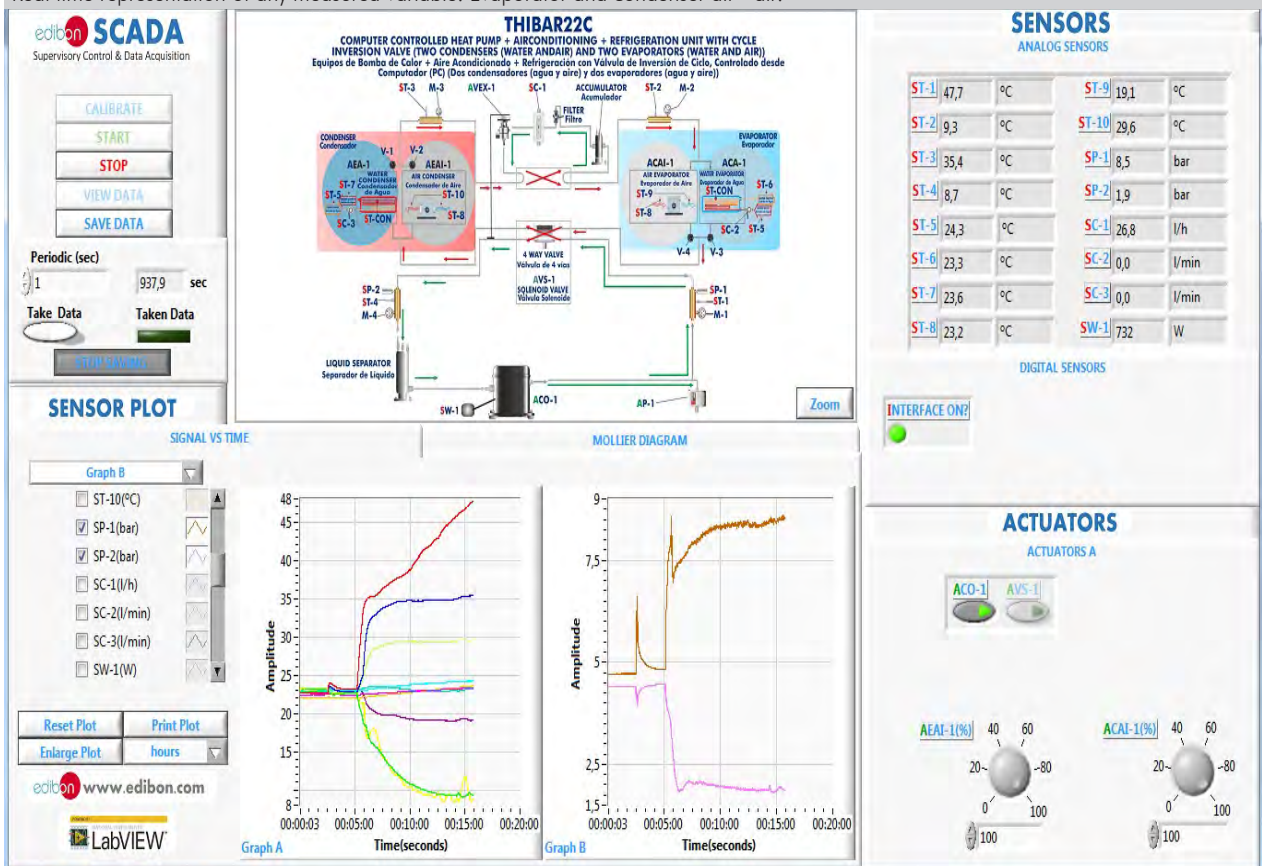

Real time representation of any measured variable. Heat pump operation. Evaporator and condenser air - water.

edibon SCADA
Supervisory Control & Data Acquisition

CALIBRATE
START
STOP
VIEW DATA
SAVE DATA

Periodic (sec)
1 / 822,8 sec

Take Data / Taken Data

SENSOR PLOT

SIGNAL VS TIME

Graph A

- ST-1(°C)
- ST-2(°C)
- ST-3(°C)
- ST-4(°C)
- ST-5(°C)
- ST-6(°C)
- ST-7(°C)

Reset Plot / Print Plot / Enlarge Plot / minutes

www.edibon.com

LabVIEW

THIBAR22C
COMPUTER CONTROLLED HEAT PUMP + AIRCONDITIONING + REFRIGERATION UNIT WITH CYCLE INVERSION VALVE (TWO CONDENSERS (WATER ANDAIR) AND TWO EVAPORATORS (WATER AND AIR))
Equipos de Bomba de Calor + Aire Acondicionado + Refrigeración con Válvula de Inversión de Ciclo, Controlado desde Computador (PC) (Dos condensadores (agua y aire) y dos evaporadores (agua y aire))

Zoom

SENSORS
ANALOG SENSORS

ST-1	38,0	°C	ST-9	27,1	°C
ST-2	6,5	°C	ST-10	30,0	°C
ST-3	35,0	°C	SP-1	8,7	bar
ST-4	5,2	°C	SP-2	1,7	bar
ST-5	18,2	°C	SC-1	25,6	l/h
ST-6	11,2	°C	SC-2	1,2	l/min
ST-7	20,0	°C	SC-3	0,0	l/min
ST-8	27,2	°C	SW-1	735	W

DIGITAL SENSORS

INTERFACE ON?

Graph A

Amplitude vs Time(seconds)

Graph B

Amplitude vs Time(seconds)

ACTUATORS
ACTUATORS A

AEAI-1(%) / ACAI-1(%)

Real time representation of any measured variable. Cooling operation. Evaporator and condenser air - water.

edibon SCADA
Supervisory Control & Data Acquisition

CALIBRATE
START
STOP
VIEW DATA
SAVE DATA

Periodic (sec)
1 / 2352,2 sec

Take Data / Taken Data

SENSOR PLOT

SIGNAL VS TIME

Graph A

- ST-1(°C)
- ST-2(°C)
- ST-3(°C)
- ST-4(°C)
- ST-5(°C)
- ST-6(°C)
- ST-7(°C)

Reset Plot / Print Plot / Enlarge Plot / minutes

www.edibon.com

LabVIEW

THIBAR22C
COMPUTER CONTROLLED HEAT PUMP + AIRCONDITIONING + REFRIGERATION UNIT WITH CYCLE INVERSION VALVE (TWO CONDENSERS (WATER ANDAIR) AND TWO EVAPORATORS (WATER AND AIR))
Equipos de Bomba de Calor + Aire Acondicionado + Refrigeración con Válvula de Inversión de Ciclo, Controlado desde Computador (PC) (Dos condensadores (agua y aire) y dos evaporadores (agua y aire))

Zoom

SENSORS
ANALOG SENSORS

ST-1	47,2	°C	ST-9	20,5	°C
ST-2	40,2	°C	ST-10	16,7	°C
ST-3	11,8	°C	SP-1	9,7	bar
ST-4	12,1	°C	SP-2	2,1	bar
ST-5	18,3	°C	SC-1	29,3	l/h
ST-6	35,3	°C	SC-2	1,3	l/min
ST-7	21,7	°C	SC-3	0,0	l/min
ST-8	21,6	°C	SW-1	771	W

DIGITAL SENSORS

INTERFACE ON?

Graph A

Amplitude vs Time(seconds)

Graph B

Amplitude vs Time(seconds)

ACTUATORS
ACTUATORS A

AEAI-1(%) / ACAI-1(%)

Real time representation of Rankine Cycle in the Mollier diagram of the coolant. Heat pump operation. Evaporator and condenser water - water.

edibon SCADA
Supervisory Control & Data Acquisition

Periodic (sec)
1 / 463,3 sec

Take Data / Taken Data

SENSOR PLOT

SIGNAL VS TIME

THIBAR22C
COMPUTER CONTROLLED HEAT PUMP + AIRCONDITIONING + REFRIGERATION UNIT WITH CYCLE INVERSION VALVE (TWO CONDENSERS (WATER ANDAIR) AND TWO EVAPORATORS (WATER AND AIR))
Equipos de Bomba de Calor + Aire Acondicionado + Refrigeración con Válvula de Inversión de Ciclo, Controlado desde Computador (PC) (Dos condensadores (agua y aire) y dos evaporadores (agua y aire))

MOLLIER DIAGRAM

SENSORS
ANALOG SENSORS

ST-1	48,9 °C	ST-9	28,0 °C
ST-2	3,8 °C	ST-10	27,9 °C
ST-3	36,1 °C	SP-1	9,7 bar
ST-4	2,6 °C	SP-2	1,4 bar
ST-5	17,4 °C	SC-1	22,8 l/h
ST-6	11,8 °C	SC-2	1,4 l/min
ST-7	30,9 °C	SC-3	1,4 l/min
ST-8	23,7 °C	SW-1	724 W

DIGITAL SENSORS

Mollier Diagram

Reset Plot / Print Plot / Enlarge Plot / minutes

www.edibon.com

LabVIEW

Real time representation of Rankine Cycle in the Mollier diagram of the coolant. Cooling operation. Evaporator and condenser water - water.

edibon SCADA
Supervisory Control & Data Acquisition

Periodic (sec)
1 / 805,2 sec

Take Data / Taken Data

SENSOR PLOT

SIGNAL VS TIME

THIBAR22C
COMPUTER CONTROLLED HEAT PUMP + AIRCONDITIONING + REFRIGERATION UNIT WITH CYCLE INVERSION VALVE (TWO CONDENSERS (WATER ANDAIR) AND TWO EVAPORATORS (WATER AND AIR))
Equipos de Bomba de Calor + Aire Acondicionado + Refrigeración con Válvula de Inversión de Ciclo, Controlado desde Computador (PC) (Dos condensadores (agua y aire) y dos evaporadores (agua y aire))

MOLLIER DIAGRAM

SENSORS
ANALOG SENSORS

ST-1	33,1 °C	ST-9	25,2 °C
ST-2	29,0 °C	ST-10	23,4 °C
ST-3	4,6 °C	SP-1	7,4 bar
ST-4	3,3 °C	SP-2	1,5 bar
ST-5	17,0 °C	SC-1	23,7 l/h
ST-6	25,4 °C	SC-2	2,4 l/min
ST-7	11,6 °C	SC-3	2,4 l/min
ST-8	23,8 °C	SW-1	713 W

DIGITAL SENSORS

Mollier Diagram


Reset Plot / Print Plot / Enlarge Plot / hours

www.edibon.com


LabVIEW

Some **real** results obtained from this Unit

Real time representation of Rankine Cycle in the Mollier diagram of the coolant.

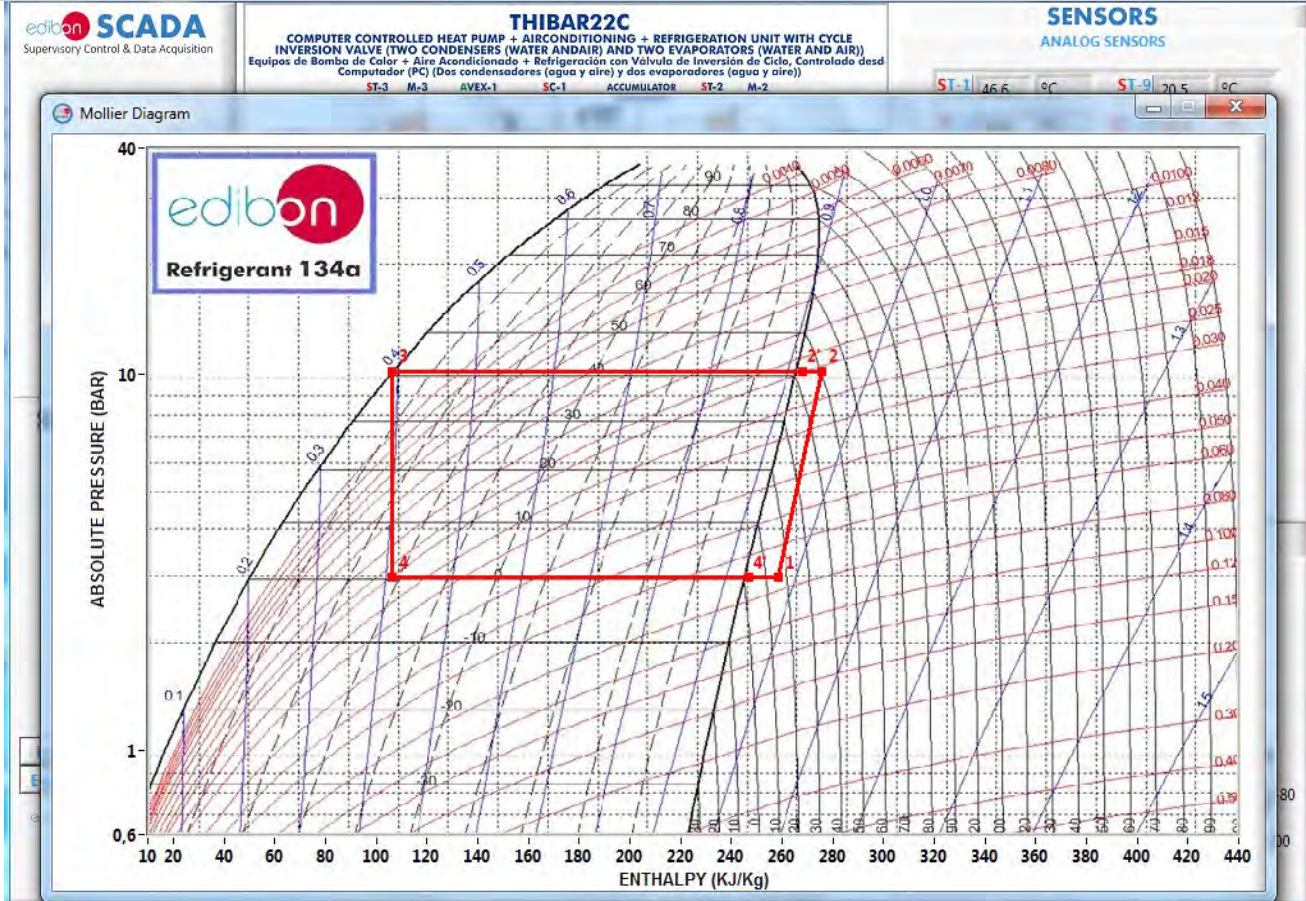

Visualization and representation of saved data.

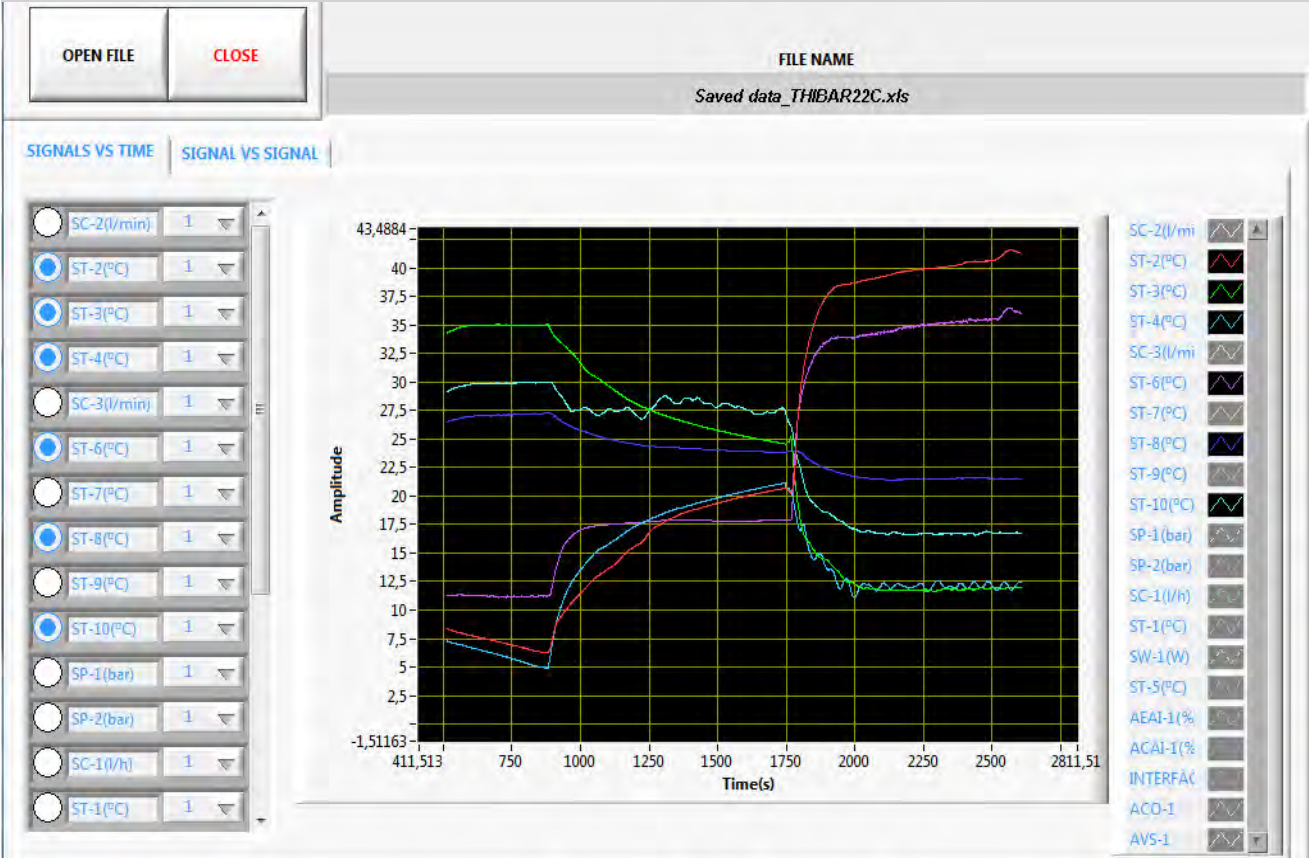

COMPLETE TECHNICAL SPECIFICATIONS (for optional items)

Additionally to the main items (1 to 6) described, we can offer, as optional, other items from 7 to 11.

All these items try to give more possibilities for:

- a) Industrial configuration. (PLC)
- b) Technical and Vocational Education configuration. (ICAI and FSS)
- c) Multipost Expansions options. (MINI ESN and ESN)

a) Industrial configuration

⑦ PLC. Industrial Control using PLC (it includes PLC-PI Module plus PLC-SOF Control Software):

-PLC-PI. PLC Module:

Metallic box.

Circuit diagram in the module front panel.

Front panel:

Digital inputs (X) and Digital outputs (Y) block:

16 Digital inputs, activated by switches and 16 LEDs for confirmation (red).

14 Digital outputs (through SCSI connector) with 14 LEDs for message (green).

Analog inputs block:

16 Analog inputs (-10 V. to + 10 V.) (through SCSI connector).

Analog outputs block:

4 Analog outputs (-10 V. to + 10 V.) (through SCSI connector).

Touch screen:

High visibility and multiple functions. Display of a highly visible status. Recipe function. Bar graph function. Flow display function. Alarm list. Multi language function. True type fonts.

Back panel:

Power supply connector. Fuse 2A. RS-232 connector to PC. USB 2.0 connector to PC.

Inside:

Power supply outputs: 24 Vdc, 12 Vdc, -12 Vdc, 12 Vdc variable.

Panasonic PLC:

High-speed scan of 0.32 μ sec. for a basic instruction.

Program capacity of 32 Ksteps, with a sufficient comment area.

Power supply input (100 to 240 V AC).

DC input: 16 (24 V DC).

Relay output: 14.

High-speed counter.

Multi-point PID control.

Digital inputs/outputs and analog inputs/outputs Panasonic modules.


Communication RS232 wire to computer (PC).

Dimensions: 490 x 330 x 310 mm. approx. (19.29 x 12.99 x 12.20 inches approx.). Weight: 30 Kg. approx. (66 pounds approx.).

-THIBAR22C/PLC-SOF. PLC Control Software:

For this particular unit, always included with PLC supply.

The software has been designed using Labview and it follows the unit operation procedure and linked with the Control Interface Box used in the Computer Controlled Reversible Heat Pump + Air Conditioning + Refrigeration with 2 Condensers and 2 Evaporators (Water/Air) (THIBAR22C).

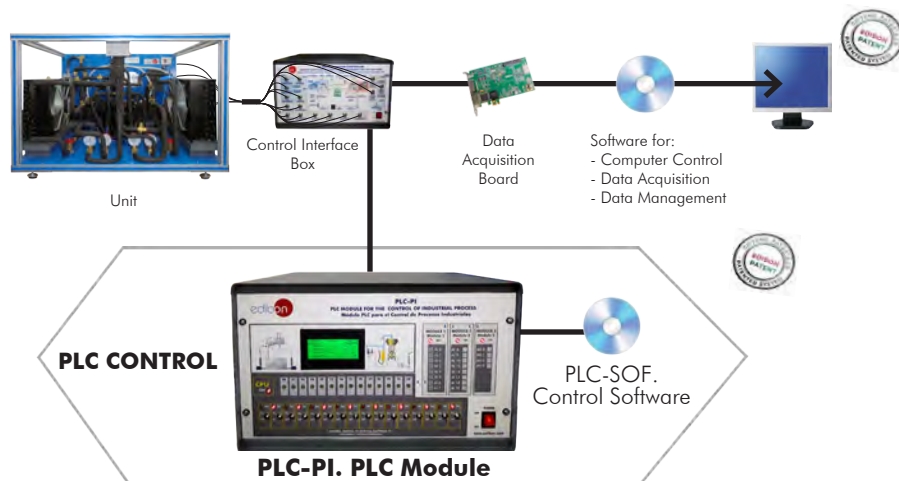

Practices to be done with PLC-PI:

- 1.- Control of the particular unit process through the control interface box without the computer.
- 2.- Visualization of all the sensors values used in the particular unit process.
- 3.- Calibration of all sensors included in the particular unit process.
- 4.- Hand on of all the actuators involved in the particular unit process.
- 5.- Realization of different experiments, in automatic way, without having in front the particular unit. (These experiments can be decided previously).
- 6.- Simulation of outside actions, in the cases do not exist hardware elements. (Example: test of complementary tanks, complementary industrial environment to the process to be studied, etc).
- 7.- PLC hardware general use.
- 8.- PLC process application for the particular unit.
- 9.- PLC structure.
- 10.- PLC inputs and outputs configuration.
- 11.- PLC configuration possibilities.
- 12.- PLC program languages.
- 13.- PLC different programming standard languages (ladder diagram (LD), structured text (ST), instructions list (IL), sequential function chart (SFC), function block diagram (FBD)).
- 14.- New configuration and development of new process.
- 15.- Hand on an established process.
- 16.- To visualize and see the results and to make comparisons with the particular unit process.
- 17.- Possibility of creating new process in relation with the particular unit.
- 18.- PLC Programming Exercises.
- 19.- Own PLC applications in accordance with teacher and student requirements.

b) Technical and Vocational Education configuration

⑧ THIBAR22C/ICAI. Interactive Computer Aided Instruction Software.

This complete software package consists of an Instructor Software (EDIBON Classroom Manager - ECM-SOF) totally integrated with the Student Software (EDIBON Student Labsoft - ESL-SOF). Both are interconnected so that the teacher knows at any moment what is the theoretical and practical knowledge of the students.

This software is optional and can be used additionally to items (1 to 6).

- ECM-SOF. EDIBON Classroom Manager (Instructor Software).

ECM-SOF is the application that allows the Instructor to register students, manage and assign tasks for workgroups, create own content to carry out Practical Exercises, choose one of the evaluation methods to check the Student knowledge and monitor the progression related to the planned tasks for individual students, workgroups, units, etc... so the teacher can know in real time the level of understanding of any student in the classroom.

Innovative features:

- User Data Base Management.
- Administration and assignment of Workgroup, Task and Training sessions.
- Creation and Integration of Practical Exercises and Multimedia Resources.
- Custom Design of Evaluation Methods.
- Creation and assignment of Formulas & Equations.
- Equation System Solver Engine.
- Updatable Contents.
- Report generation, User Progression Monitoring and Statistics.

- ESL-SOF. EDIBON Student Labsoft (Student Software).

ESL-SOF is the application addressed to the Students that helps them to understand theoretical concepts by means of practical exercises and to prove their knowledge and progression by performing tests and calculations in addition to Multimedia Resources. Default planned tasks and an Open workgroup are provided by EDIBON to allow the students start working from the first session. Reports and statistics are available to know their progression at any time, as well as explanations for every exercise to reinforce the theoretically acquired technical knowledge.

Innovative features:

- Student Log-In & Self-Registration.
- Existing Tasks checking & Monitoring.
- Default contents & scheduled tasks available to be used from the first session.
- Practical Exercises accomplishment by following the Manual provided by EDIBON.
- Evaluation Methods to prove your knowledge and progression.
- Test self-correction.
- Calculations computing and plotting.
- Equation System Solver Engine.
- User Monitoring Learning & Printable Reports.
- Multimedia-Supported auxiliary resources.

For more information see ICAI catalogue. Click on the following link:
www.edibon.com/en/files/expansion/ICAI/catalog

Instructor Software



ECM-SOF. EDIBON Classroom Manager (Instructor Software) Application Main Screen


ECAL. EDIBON Calculations Program Package - Formula Editor Screen



ERS. EDIBON Results & Statistics Program Package - Student Scores Histogram


ETTE. EDIBON Training Test & Exam Program Package - Main Screen with Numeric Result Question

Student Software

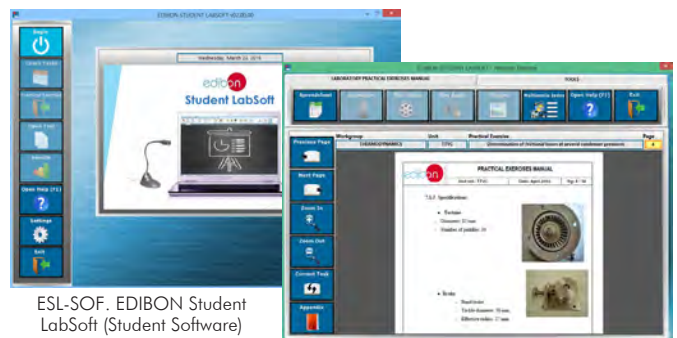

ESL-SOF. EDIBON Student LabSoft (Student Software) Application Main Screen

EPE. EDIBON Practical Exercise Program Package Main Screen

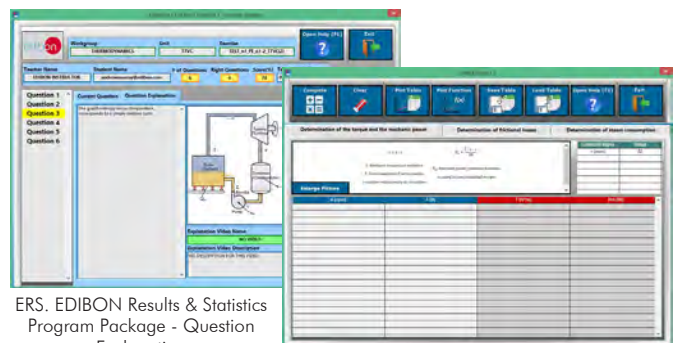

ERS. EDIBON Results & Statistics Program Package - Question Explanation

ECAL. EDIBON Calculations Program Package Main Screen

⑨ **THIBAR22C/FSS. Faults Simulation System.**

Faults Simulation System (FSS) is a Software package that simulates several faults in any EDIBON Computer Controlled Unit. It is useful for Technical and Vocational level.


The "FAULTS" mode consists in causing several faults in the unit normal operation. The student must find them and solve them. There are several kinds of faults that can be grouped in the following sections:

- Faults affecting the sensors measurement:
 - An incorrect calibration is applied to them.
 - Non-linearity.
- Faults affecting the actuators:
 - Actuators channels interchange at any time during the program execution.
 - Response reduction of an actuator.
- Faults in the controls execution:
 - Inversion of the performance in ON/OFF controls.
 - Reduction or increase of the calculated total response.
 - The action of some controls is annulled.
- On/off faults:
 - Several on/off faults can be included.

For more information see FSS catalogue. Click on the following link:

www.edibon.com/en/files/expansion/FSS/catalog

Example of some screens

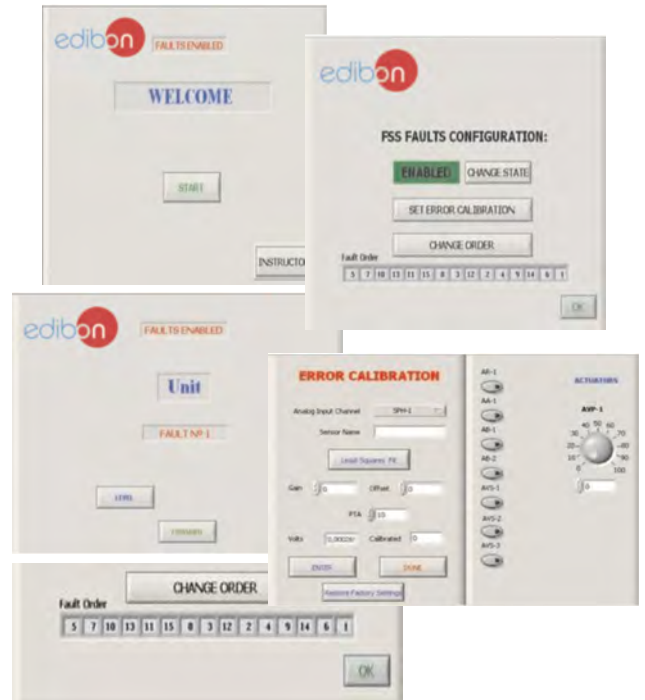

c) Multipost Expansions options

⑩ **MINI ESN. EDIBON Mini Scada-Net System for being used with EDIBON Teaching Units.**

MINI ESN. EDIBON Mini Scada-Net System allows up to 30 students to work with a Teaching Unit in any laboratory, simultaneously. It is useful for both, Higher Education and/or Technical and Vocational Education.

The MINI ESN system consists of the adaptation of any EDIBON Computer Controlled Unit with SCADA integrated in a local network.

This system allows to view/control the unit remotely, from any computer integrated in the local net (in the classroom), through the main computer connected to the unit. Then, the number of possible users who can work with the same unit is higher than in an usual way of working (usually only one).

Main characteristics:

- It allows up to 30 students to work simultaneously with the EDIBON Computer Controlled Unit with SCADA, connected in a local net.
- Open Control + Multicontrol + Real Time Control + Multi Student Post.
- Instructor controls and explains to all students at the same time.
- Any user/student can work doing "real time" control/multicontrol and visualisation.
- Instructor can see in the computer what any user/student is doing in the unit.
- Continuous communication between the instructor and all the users/students connected.

Main advantages:

- It allows an easier and quicker understanding.
- This system allows you can save time and cost.
- Future expansions with more EDIBON Units.

For more information see MINI ESN catalogue. Click on the following link:

www.edibon.com/en/files/expansion/MINI-ESN/catalog


⑪ **ESN. EDIBON Scada-Net Systems.**

This unit can be integrated, in the future, into a Complete Laboratory with many Units and many Students.

For more information see ESN catalogue. Click on the following link:

www.edibon.com/en/files/expansion/ESN/catalog


MINI ESN.
EDIBON Mini Scada-Net System



1 UNIT= up to 30 STUDENTS can work simultaneously

Computer Controlled Reversible Heat Pump + Air Conditioning + Refrigeration with 2 Condensers and 2 Evaporators (Water/Air) (THIBAR22C)


Control Interface Box




Software for:

- Computer Control
- Data Acquisition
- Data Management


Instructors' Central Computer



Mini Scada-Net Software



LOCAL NET FOR 30 STUDENTS

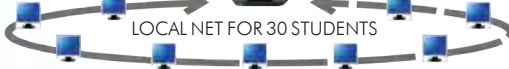

OPEN CONTROL + MULTICONTROL + REAL TIME CONTROL + MULTI STUDENT POST

Note: The MINI ESN system can be used with any EDIBON computer controlled unit

ORDER INFORMATION

Main items (always included in the supply)

Minimum supply always includes:

- ① **Unit: THIBAR22C. Computer Controlled Reversible Heat Pump + Air Conditioning + Refrigeration with 2 Condensers and 2 Evaporators (Water/Air).**
- ② **THIBAR22C/CIB. Control Interface Box.**
- ③ **DAB. Data Acquisition Board.**
- ④ **THIBAR22C/CCSOF. Computer Control + Data Acquisition + Data Management Software.**
- ⑤ **Cables and Accessories**, for normal operation.
- ⑥ **Manuals.**

***IMPORTANT:** Under THIBAR22C we always supply all the elements for immediate running as 1, 2, 3, 4, 5 and 6.

Optional items (supplied under specific order)

a) Industrial configuration

- ⑦ PLC. Industrial Control using PLC (it includes PLC-PI Module plus PLC-SOF Control Software):
 - PCL-PI. PLC Module.
 - THIBAR22C/PLC-SOF. PLC Control Software.

b) Technical and Vocational Education configuration

- ⑧ THIBAR22C/ICAL. Interactive Computer Aided Instruction Software.
- ⑨ THIBAR22C/FSS. Faults Simulation System.

c) Multipost Expansions options

- ⑩ MINI ESN. EDIBON Mini Scada-Net System for being used with EDIBON Teaching Units.
- ⑪ ESN. EDIBON Scada-Net Systems.

① THIBAR22C. Unit:

Bench-top unit.
 Anodized aluminum frame and panels made of painted steel.
 Main metallic elements made of stainless steel.
 Diagram in the front panel with distribution of the elements similar to the real one.
 Hermetic compressor, power: 3/8 HP, computer controlled.
 Water condenser, capacity: 1580 W.
 Air condenser, computer controlled.
 High pressure control.
 Storage vessel to prevent a wrong operation of the compressor, capacity: 0.5 l.
 Filter for the working fluid.
 Expansion valve.
 Water evaporator, capacity: 1580 W.
 Air evaporator, computer controlled.
 Liquid separator to prevent liquid from entering the compressor, capacity: 0.73 l.
 Four manometers, situated at important points of the unit.
 Ten temperature sensors (four sensors measure the cooling temperature, three sensors measure the water temperature, three sensors measure the air temperature):
 "J" type temperature sensor (compressor outlet).
 "J" type temperature sensor (condenser outlet/evaporator inlet).
 "J" type temperature sensor (evaporator inlet/condenser outlet).
 "J" type temperature sensor (compressor inlet).
 "J" type temperature sensor (water inlet).
 "J" type temperature sensor (condenser water outlet).
 "J" type temperature sensor (evaporator water outlet).
 "J" type temperature sensor (air inlet).
 "J" type temperature sensor (condenser air outlet).
 "J" type temperature sensor (evaporator air outlet).
 Three flow sensors:
 Cooling flow sensor, range: 5 – 60 l/h.
 Water flow sensor (water condenser), range: 0.25 – 6.5 l/min.
 Water flow sensor (water evaporator), range: 0.25 – 6.5 l/min.
 Two pressure sensors:
 Cooling pressure sensor (compressor outlet), range: 0 – 25 bar.
 Cooling pressure sensor (compressor inlet), range: 0 – 10 bar.
 Compressor wattmeter.
 Four-way valve (solenoid valve) for reversibility the cycle.
 Four valves to direct the coolant through the condensers and evaporators.
 Enthalpy diagram of the R134a coolant.
 The complete unit includes as well:
 Advanced Real-Time SCADA.
 Open Control + Multicontrol + Real-Time Control.
 Specialized EDIBON Control Software based on LabVIEW.
 National Instruments Data Acquisition board (250 KS/s, kilo samples per second).
 Calibration exercises, which are included, teach the user how to calibrate a sensor and the importance of checking the accuracy of the sensors before taking measurements.
 Projector and/or electronic whiteboard compatibility allows the unit to be explained and demonstrated to an entire class at one time.
 Capable of doing applied research, real industrial simulation, training courses, etc.
 Remote operation and control by the user and remote control for EDIBON technical support, are always included.
 Totally safe, utilizing 4 safety systems (Mechanical, Electrical, Electronic & Software).
 Designed and manufactured under several quality standards.
 Optional ICAI software to create, edit and carry out practical exercises, tests, exams, calculations, etc. Apart from monitoring user's knowledge and progress reached.
 This unit has been designed for future expansion and integration. A common expansion is the EDIBON Scada-Net (ESN) System which enables multiple students to simultaneously operate many units in a network.

② THIBAR22C/CIB. Control Interface Box:

The Control Interface Box is part of the SCADA system.
 Control interface box with process diagram in the front panel.
 The unit control elements are permanently computer controlled.
 Simultaneous visualization in the computer of all parameters involved in the process.
 Calibration of all sensors involved in the process.
 Real time curves representation about system responses.
 All the actuators' values can be changed at any time from the keyboard allowing the analysis about curves and responses of the whole process.
 Shield and filtered signals to avoid external interferences.
 Real time computer control with flexibility of modifications from the computer keyboard of the parameters, at any moment during the process.
 Real time computer control for parameters involved in the process simultaneously.
 Open control allowing modifications, at any moment and in real time, of parameters involved in the process simultaneously.
 Three safety levels, one mechanical in the unit, another electronic in the control interface and the third one in the control software.

③ DAB. Data Acquisition Board:

The Data Acquisition board is part of the SCADA system.
 PCI Express Data acquisition board (National Instruments) to be placed in a computer slot.
 Analog input: Channels= 16 single-ended or 8 differential. Resolution= 16 bits, 1 in 65536. Sampling rate up to: 250 KS/s (kilo samples per second).
 Analog output: Channels=2. Resolution=16 bits, 1 in 65536.
 Digital Input/Output: Channels=24 inputs/outputs.
 The Data Acquisition board model may change at any moment, providing the same or better features than those required for the unit.

④ THIBAR22C/CCSOF. Computer Control + Data Acquisition+ Data Management Software:

The three softwares are part of the SCADA system.
 Compatible with the industry standards.
 Flexible, open and multicontrol software, developed with actual windows graphic systems, acting simultaneously on all process parameters.
 Management, processing, comparison and storage of data.
 Sampling velocity up to 250 KS/s (kilo samples per second).
 Calibration system for the sensors involved in the process.
 It allows the registration of the alarms state and the graphic representation in real time.
 Open software, allowing the teacher to modify texts, instructions. Teacher's and student's passwords to facilitate the teacher's control on the student, and allowing the access to different work levels.
 This unit allows the 30 students of the classroom to visualize simultaneously all the results and the manipulation of the unit, during the process, by using a projector or an electronic whiteboard.

⑤ Cables and Accessories, for normal operation.**⑥ Manuals:**

This unit is supplied with 8 manuals: Required Services, Assembly and Installation, Interface and Control Software, Starting-up, Safety, Maintenance, Calibration & Practices Manuals.

Exercises and Practical Possibilities to be done with the Main Items

- 1.- Determination of COP (coefficient of performance) of a heat pump. Water as heat source. (Water-water heat pump).
 - 2.- Determination of COP (coefficient of performance) of a heat pump. Air as heat source. (Water-air heat pump).
 - 3.- Determination of COP (coefficient of performance) of a heat pump. Air as heat source. (Air-air heat pump).
 - 4.- Determination of COP (coefficient of performance) of a heat pump. Water as heat source. (Air-water heat pump).
 - 5.- Representation of performance curves of the heat pump with different inlet and outlet temperatures. Water as heat source. (Water-water heat pump).
 - 6.- Representation of performance curves of the heat pump with different inlet and outlet temperatures. Air as heat source. (Water-air heat pump).
 - 7.- Representation of performance curves of the heat pump with different inlet and outlet temperatures. Water as heat source. (Air-water heat pump).
 - 8.- Representation of performance curves of the heat pump with different inlet and outlet temperatures. Air as heat source. (Air-air heat pump).
 - 9.- Lay out of the refrigeration compression cycle in a diagram P-H and comparison with the ideal cycle. Water as heat source. (Water-water heat pump).
 - 10.- Lay out of the refrigeration compression cycle in a diagram P-H and comparison with the ideal cycle. Air as heat source. (Water-air heat pump).
 - 11.- Lay out of the refrigeration compression cycle in a diagram P-H and comparison with the ideal cycle. Water as heat source. (Air-water heat pump).
 - 12.- Lay out of the refrigeration compression cycle in a diagram P-H and comparison with the ideal cycle. Air as heat source. (Air-air heat pump).
 - 13.- Representation of the performance curves of the heat pump based on the properties of the refrigerant and at different condensation and evaporation temperatures. Water as heat source. (Water-water heat pump).
 - 14.- Representation of the performance curves of the heat pump based on the properties of the refrigerant and at different condensation and evaporation temperatures. Air as heat source. (Water-air heat pump).
 - 15.- Representation of the performance curves of the heat pump based on the properties of the refrigerant and at different condensation and evaporation temperatures. Water as heat source. (Air-water heat pump).
 - 16.- Representation of the performance curves of the heat pump based on the properties of the refrigerant and at different condensation and evaporation temperatures. Air as heat source. (Air-air heat pump).
 - 17.- Practices with cycle inversion.
- Additional practical possibilities:
- 18.- Sensors calibration.
- Other possibilities to be done with this Unit:
- 19.- Many students view results simultaneously.
To view all results in real time in the classroom by means of a projector or an electronic whiteboard.
 - 20.- Open Control, Multicontrol and Real Time Control.
This unit allows intrinsically and/or extrinsically to change the span, gains; proportional, integral, derivative parameters; etc, in real time.
 - 21.- The Computer Control System with SCADA allows a real industrial simulation.
 - 22.- This unit is totally safe as uses mechanical, electrical/electronic, and software safety devices.
 - 23.- This unit can be used for doing applied research.
 - 24.- This unit can be used for giving training courses to Industries even to other Technical Education Institutions.
 - 25.- Control of the THIBAR22C unit process through the control interface box without the computer.
 - 26.- Visualization of all the sensors values used in the THIBAR22C unit process.
- By using PLC-PI additional 19 more exercises can be done.
 - Several other exercises can be done and designed by the user.

TENDER SPECIFICATIONS (for optional items)

a) Industrial configuration

⑦ **PLC. Industrial Control using PLC** (it includes PLC-PI Module plus PLC-SOF Control Software):

-PLC-PI. PLC Module:

Metallic box.

Circuit diagram in the module front panel.

Digital inputs (X) and Digital outputs (Y) block: 16 Digital inputs. 14 Digital outputs.

Analog inputs block: 16 Analog inputs.

Analog outputs block: 4 Analog outputs.

Touch screen.

Panasonic PLC:

High-speed scan of 0.32 μ sec. Program capacity of 32 Ksteps. High-speed counter. Multi-point PID control.

Digital inputs/outputs and analog inputs/outputs Panasonic modules.

-THIBAR22C/PLC-SOF. PLC Control Software:

For this particular unit, always included with PLC supply.

Practices to be done with PLC-PI:

- 1.- Control of the particular unit process through the control interface box without the computer.
- 2.- Visualization of all the sensors values used in the particular unit process.
- 3.- Calibration of all sensors included in the particular unit process.
- 4.- Hand on of all the actuators involved in the particular unit process.
- 5.- Realization of different experiments, in automatic way, without having in front the particular unit. (These experiments can be decided previously).
- 6.- Simulation of outside actions, in the cases do not exist hardware elements. (Example: test of complementary tanks, complementary industrial environment to the process to be studied, etc).
- 7.- PLC hardware general use.
- 8.- PLC process application for the particular unit.
- 9.- PLC structure.
- 10.- PLC inputs and outputs configuration.
- 11.- PLC configuration possibilities.
- 12.- PLC program languages.
- 13.- PLC different programming standard languages (ladder diagram (LD), structured text (ST), instructions list (IL), sequential function chart (SFC), function block diagram (FBD)).
- 14.- New configuration and development of new process.
- 15.- Hand on an established process.
- 16.- To visualize and see the results and to make comparisons with the particular unit process.
- 17.- Possibility of creating new process in relation with the particular unit.
- 18.- PLC Programming Exercises.
- 19.- Own PLC applications in accordance with teacher and student requirements.

b) Technical and Vocational Education configuration

⑧ **THIBAR22C/ICAI. Interactive Computer Aided Instruction Software.**

This complete software package consists of an Instructor Software (EDIBON Classroom Manager - ECM-SOF) totally integrated with the Student Software (EDIBON Student Labsoft - ESL-SOF). Both are interconnected so that the teacher knows at any moment what is the theoretical and practical knowledge of the students.

- ECM-SOF. EDIBON Classroom Manager (Instructor Software).

ECM-SOF is the application that allows the Instructor to register students, manage and assign tasks for workgroups, create own content to carry out Practical Exercises, choose one of the evaluation methods to check the Student knowledge and monitor the progression related to the planned tasks for individual students, workgroups, units, etc...so the teacher can know in real time the level of understanding of any student in the classroom.

Innovative features:

- User Data Base Management.
- Administration and assignment of Workgroup, Task and Training sessions.
- Creation and Integration of Practical Exercises and Multimedia Resources.
- Custom Design of Evaluation Methods.
- Creation and assignment of Formulas & Equations.
- Equation System Solver Engine.
- Updatable Contents.
- Report generation, User Progression Monitoring and Statistics.

- ESL-SOF. EDIBON Student Labsoft (Student Software).

ESL-SOF is the application addressed to the Students that helps them to understand theoretical concepts by means of practical exercises and to prove their knowledge and progression by performing tests and calculations in addition to Multimedia Resources. Default planned tasks and an Open workgroup are provided by EDIBON to allow the students start working from the first session. Reports and statistics are available to know their progression at any time, as well as explanations for every exercise to reinforce the theoretically acquired technical knowledge.

Innovative features:

- Student Log-In & Self-Registration.
- Existing Tasks checking & Monitoring.
- Default contents & scheduled tasks available to be used from the first session.
- Practical Exercises accomplishment by following the Manual provided by EDIBON.
- Evaluation Methods to prove your knowledge and progression.
- Test self-correction.
- Calculations computing and plotting.
- Equation System Solver Engine.
- User Monitoring Learning & Printable Reports.
- Multimedia-Supported auxiliary resources.

⑨ **THIBAR22C/FSS. Faults Simulation System.**

Faults Simulation System (FSS) is a Software package that simulates several faults in any EDIBON Computer Controlled Unit. The "FAULTS" mode consists in causing several faults in the unit normal operation. The student must find them and solve them. There are several kinds of faults that can be grouped in the following sections:

Faults affecting the sensors measurement:

- An incorrect calibration is applied to them.
- Non-linearity.

Faults affecting the actuators:

- Actuators channels interchange at any time during the program execution.
- Response reduction of an actuator.

Faults in the controls execution:

- Inversion of the performance in ON/OFF controls.
- Reduction or increase of the calculated total response.
- The action of some controls is annulled.

On/off faults:

- Several on/off faults can be included.

c) Multipost Expansions options

⑩ **MINI ESN. EDIBON Mini Scada-Net System for being used with EDIBON Teaching Units.**

MINI ESN. EDIBON Mini Scada-Net System allows up to 30 students to work with a Teaching Unit in any laboratory, simultaneously.

The MINI ESN system consists of the adaptation of any EDIBON Computer Controlled Unit with SCADA integrated in a local network.

This system allows to view/control the unit remotely, from any computer integrated in the local net (in the classroom), through the main computer connected to the unit.

Main characteristics:

- It allows up to 30 students to work simultaneously with the EDIBON Computer Controlled Unit with SCADA, connected in a local net.
- Open Control + Multicontrol + Real Time Control + Multi Student Post.
- Instructor controls and explains to all students at the same time.
- Any user/student can work doing "real time" control/multicontrol and visualisation.
- Instructor can see in the computer what any user/student is doing in the unit.
- Continuous communication between the instructor and all the users/students connected.

Main advantages:

- It allows an easier and quicker understanding.
- This system allows you can save time and cost.
- Future expansions with more EDIBON Units.

The system basically will consist of:

This system is used with a Computer Controlled Unit.

- Instructor's computer.
- Students' computers.
- Local Network.
- Unit-Control Interface adaptation.
- Unit Software adaptation.
- Webcam.
- MINI ESN Software to control the whole system.
- Cables and accessories required for a normal operation.

* Specifications subject to change without previous notice, due to the convenience of improvement of the product.



C/ Julio Cervera, 10-12-14. Móstoles Tecnológico.
28935 MÓSTOLES. (Madrid). ESPAÑA - SPAIN.
Tel.: 34-91-6199363 Fax: 34-91-6198647

E-mail: edibon@edibon.com Web: www.edibon.com

Edition: ED01/20
Date: September/2020

REPRESENTATIVE:

